

Critical Environment Solutions

HEPA & ULPA Filters Handling Instructions



DO NOT use a knife to open the box or filter bag.



DO NOT Stack filters or place anything on the filter media.



DO NOT handle or manipulate the filter media or gel seal.



DO Lift filters only vertically and by the corners. Use two people for filters 2'x3' or larger.

HANDLING INSTRUCTIONS

- Keep filters boxed and bagged.
- Lift and carry the filter vertically. Handle with clean, dry hands. Hold filters by the corners of the frame.
- Use two people to lift filters larger than 2'x2' to avoid twisting the frame.
- DO NOT use a knife to open filter boxes and bags.
- DO NOT touch filter gaskets, gel seals, or the filter media.
- DO NOT twist the filter frame.
- DO NOT drop or knock the filter.

STORAGE INSTRUCTIONS

- Keep filters in an upright vertical orientation.
- DO NOT get filters wet.
- DO NOT stack filters or store anything on top of filters or filter boxes.
- DO NOT store filters in direct sunlight

INSTALLATION INSTRUCTIONS

- Replace the filter if your maintenance program calls for it, filter integrity is compromised, face velocity falls below acceptable levels, pressure drop across the filter has doubled since install, or the filter has reached 5 years of use or 10 years of age.
- Wear personal protective equipment when installing filters; Goggles, masks, gloves, and coveralls.
- Filters are not washable or reusable. Spent filters should be discarded in accordance with local regulation.