


Softwall Modular Cleanroom

© Copyright 2021 Terra Universal Inc. All rights reserved. • Revised Sept. 2022

Your Comprehensive Equipment Source

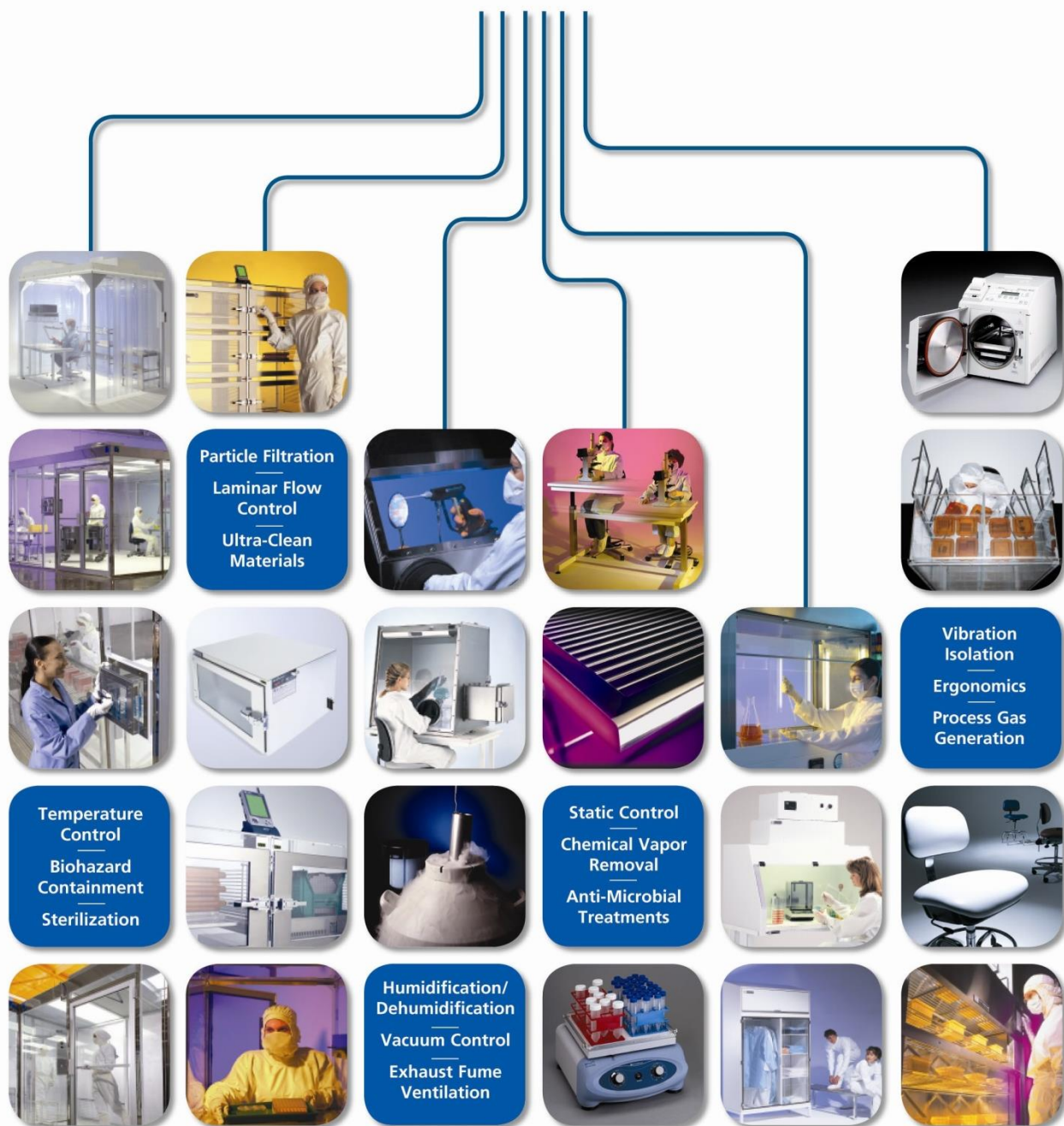



Softwall Modular Cleanroom

Table of Contents

Section 1.0 – Description

Description	3
Frame	3
Fan Filter Units.....	4
Light Panel.....	5
Blank Ceiling Panels	5
Softwall Panels & Strip Shields	5
Casters	6

Section 2.0 –Installation

Site Preparation	6
Required Installation Equipment	6
Components List	6
Cleanroom Frame Components.....	8
Ceiling Components.....	8
Support Bar Components.....	9
Top Trimming Components.....	9
Curtain Panel Components	10
Frame Assembly	10
Seismic Anchoring	13
Joists and Divider Installation.....	14
Use of the Spacer	17
Support Bar Installation (For Cleanrooms from 10 ft. to 12 ft.).....	19
Installing 2 Support Bars (For Rooms Bigger Than 12 ft.).....	21
Top Trimming Installation.....	22
Installing the Ceiling Modules	24
Curtain Panel Installation	26

Section 3.0 – Electrical Wiring

Installing the PDM.....	28
Operation Status Indicator Lights Wiring.....	31
Fan Filter Unit Hard Wiring.....	32
Fan Filter Unit Speed Configuration with PDM.....	33

Section 4.0 – Operation & Maintenance

Filter Replacement.....	35
Cleanroom Maintenance	37

Section 5.0 – Warranty

Warranty Information	38
----------------------------	----

Section 6.0 – Spare Parts & Accessories

Spare Parts & Accessories	39
---------------------------------	----



Softwall Modular Cleanroom

Proprietary Notice

This manual pertains to proprietary devices manufactured by Terra Universal, Inc. Neither this document nor any portion of it may be reproduced in any way without prior written permission from Terra Universal.

Terra Universal makes no warranties applying to information contained in this manual or its suitability for any implied or inferred purpose. Terra Universal shall not be held liable for any errors this manual contains or for any damages that result from its use.

Safety Notice

A thorough familiarity with all operating guidelines is essential to safe operation of the product. Failure to observe safety precautions could result in poor performance, damage to the system or other property, or serious bodily injury or death.

The following symbols are intended to call your attention to two levels of hazard involved in operation:



Cautions are used when failure to observe instructions could result in significant damage to equipment.



Warnings are used when failure to observe instructions or precautions could result in injury or death.

The information presented here is subject to change without notice.



IMPORTANT SAFETY NOTICE

Terra Universal cleanrooms are not designed to support more weight than the blower modules and lighting fixtures originally installed. Ceiling grid beams are no load-bearing and will not support personnel or other additional loads. Placing added weight on the ceiling grid may result in serious damage to the cleanroom and its occupants.

Safety notices supplied by Terra Universal must be affixed prominently on each side of the cleanroom grid.

1.0 Description

Terra Universal's Softwall Modular Cleanrooms provide an economical, versatile solution to clean manufacturing requirements in many industries. Portable, expandable, and easy to assemble/disassemble, this system features ceiling HEPA fan filter units that direct a continuous flow of particle-free air through the enclosure. It is completely free standing and requires no external bracing.

Frame

This system consists of a steel support frame, ceiling mounted filtration and illumination system, flexible side panels and a variety of other access options. The softwall cleanroom support frame is fabricated of interlocking 2.5" powder-coated square steel or stainless steel uprights, cross sections, and ceiling members.



Ceiling members are connected to steel T-bars, forming the 2' X 4' bays that support filters and lights. Corner reinforcement gussets ensure a rigid structure. Optional heavy-duty casters allow easy repositioning after installation. Standard height is seven feet measured from filter face to floor (8' 3" overall), allowing installation below a standard nine foot ceiling. Custom ceiling heights up to ten feet are available. System includes a control panel for operation of fan filter units, lights, and other selected options.

Type	Cat. #
Powder-Coated Steel	6600-60
304 Stainless Steel	6600-61



Softwall Modular Cleanroom

Fan Filter Units

Fan Filter Units direct a vertical laminar flow of filtered air downward through the enclosed cleanroom area.

Each includes a 650 CFM impeller blower (rated at 90 FPM with filter load) mounted in an aluminum housing with a plenum design that ensures ample air velocity across the entire face of the filter. A HEPA (high efficiency particulate air) filter installed inside the housing is rated 99.99% efficient at 0.3um particles. The filtration medium consists of micro porous polyurethane mini-pleats held in place by strong, rigid plastic separators that keep the medium from nesting. This design channels airflow with optimal efficiency to reduce resistance. The filter is sealed into the sturdy powder-coated steel frame with a fire-retardant, non-outgassing adhesive. On an optional basis, an ULPA (ultra-low penetration air) filter, rated 99.9995% efficient at 0.12um particles, may be substituted for the HEPA filter. Power to the Fan Filter Units is controlled by a master ON/OFF switch located on the cleanroom control panel. All 120VAC units and 220VAC, 60Hz units are UL listed. CE-marked models are available for 220VAC, 50Hz operations.

Filter Fan Units (HP: high speed; MS: medium speed; LS: low speed)										
Model	Dimensions Inches (mm)	Wt. Lbs. (Kg)	Avg. CFM @ MS	Airflow ft/min (m/s)			Run Amps Watts @ 60Hz			Power
				HS	MS	LS	HS	MS	LS	
2 x 4	23.63 x 47.63 x 12.51 (600 x 1210 x 318)	71 (32)	717 (1218)	115 (.58)	102 (.51)	93 (.47)	4.3A (512W)	3.5A (416W)	3.3A (393W)	120VAC, 60 Hz 220VAC, 50/60 Hz
2 x 3	23.63 x 35.63 x 12.51 (600 x 905 x 318)	53 (24)	602 (1023)	121 (.61)	116 (.59)	106 (.54)	4.1A (500W)	3.1A (378W)	2.7A (329W)	120VAC, 60 Hz 220VAC, 50/60 Hz
2 x 2	23.63 x 23.63 x 12.51 (600 x 600 x 318)	44 (20)	558 (948)	172 (.87)	166 (.84)	162 (.82)	3.9W (472W)	2.8A (339W)	2.4A (290W)	120VAC, 60 Hz 220VAC, 50/60 Hz

Model: Terra Universal WhisperFlow Fan Filter Unit



If your order included UL documentation (Cat. # 6600-33), your cleanroom will include labels indicating the use of UL-listed components on one of the triangular gussets (see photo to the right).

Recommended Fan Filter Configurations (based on ceiling height of 8 feet or less)

Recommended Fan Filter Configurations							
Nominal Dimensions	No. of Ceiling Bays	No. of Filter Modules to meet Fed Std 209 Class Standards					
		1	10	100	1,000	10,000	100,000
8' x 8'	8	5-8	4-8	3-6	2-3	1-2	1
8' x 12'	12	7-12	6-12	4-8	3-5	2-3	1-2
8' x 16'	16	10-16	8-16	6-11	4-6	2-4	1-3
12' x 12'	18	11-18	9-18	6-13	5-7	3-5	1-3
12' x 16'	24	14-24	12-24	8-17	6-10	4-6	2-4
16' x 16'	32	19-32	16-32	11-22	8-13	5-8	2-5
20' x 20'	50	30-50	25-50	18-35	13-20	8-13	3-8





Softwall Modular Cleanroom

Light Panels



Each light panel is externally mounted to ensure effective illumination of the work area without interfering with the controlled air stream. Each module is controlled by a master switch mounted on the main Cleanroom Control Panel.

Cat. # 3800-41B

Power Requirements

USA: 115VAC/60Hz, 1.50A Each

International: 277VAC/50Hz, 0.65A Each

Blank Ceiling Panels

All remaining ceiling grids are covered with blank panels, which can be removed to allow installation of additional fan filter units or illuminator modules. Panels are made of white polypropylene, acrylic or steel depending on customer selection. All ceiling modules rest against the ceiling grid frame to form a tight seal along the perimeters.

Softwall Panels & Strip Shields

Softwall panels are fabricated of heavy-duty, blur-free 40 mil, vinyl or polyurethane—double the industry standard—ensuring enhanced tear resistance and eliminating the need for unsightly sewn sash weights. Strip shields, made of 60-mil material, come standard in 8" widths, with 2" overlap, for easy entrance or exiting. Polyurethane shields are 12" W, 40 mil thick.

Panels and strips are available in three materials, in either clear or frosted appearance.

- Cleanroom-Grade Vinyl is fire retardant and low outgassing
- Static-Dissipative Vinyl is manufactured with permanent built in static inhibitors that unlike topical treatments, will not wear off.
- Non-outgassing panels are made from inherently static-dissipative polyurethane that provides permanent, humidity-independent ESD protection for sensitive manufacturing conditions. Polyurethane contains no plasticizer, making it virtually non-outgassing.



Note: Softwall panels are notched to ensure a snug fit at all corners. In standard installations, panels are supplied in lengths of four feet. Where two panels meet, the panels overlap by 6". To prevent personnel from parting panels at these seams, order optional panel seaming, Dual-Lock Velcro, magnetic closures or attachment clips below.

Panel Options	Cat. #
Dual – Lock Velcro Strips	6704-05
Flexible Magnet Strip Self-Closures	6704-06
Snap – Tite™ Clips	6704-07
Curtain Seaming	6600-65



Softwall Modular Cleanroom

Casters



Stainless steel swivel casters feature nonshedding polyurethane wheels. Capacity: 300 lbs. each. Order one per vertical post. Casters raise the frame 7.125". If casters are ordered, the frame will be made shorter to keep the floor to filter face dimension to specification. Order one per vertical post.

Cat. # 6704-08



Caution: Casters are intended to allow movement only on smooth floors and require that personnel support each cleanroom upright during motion.

2.0 Installation

Unpack all system components and check for damaged. Any damage should be reported to the shipping company immediately.

Site Preparation

- A. Site preparation requirements should be observed for any installation performed by Terra Universal technicians.
- B. Customer facility area where the cleanroom is to be installed must afford a minimum clearance of one foot (including fixtures ducts and pipes) on all sides and at least two feet clearance between FFU inlet and ceiling.
- C. Customers must provide permanent electrical connection from supply panel to Terra Cleanroom electrical junction boxes (one power line per junction box) in conformance with local electrical code, as well as any vacuum, air, H₂O, sprinkler, or nitrogen connections required for the cleanroom.
- D. Customers must give advance notice of dates and times for Terra personnel to perform the installation. Short notice may result in higher fees for travel and accommodations.
- E. Customers are to provide utilities, installation power and removal of any packing material.
- F. Because installation requires unpacking and assembling components, customers are to ensure an adequate staging area for parts and equipment adjacent to the assembly area, clear and ready for work.
- G. Customers are to inform Terra Universal in advance of any requirements for security, escorts, special training, badges, work hours, parking areas or special identification and how to obtain all such required permissions and related items.

Required Installation Equipment

- A. Handling cleanroom shipping crates, which generally measure over 300" long and weigh well over 1,000 pounds, requires at least one forklift. If crates must be moved through narrow aisles or entrances, two forklifts or pallet jacks are recommended, one to support each end.
- B. Unloading crates from the truck is much easier if you have a truck-high loading dock. Without such a dock, you will need at least one forklift and a support to brace one end while the forklift is positioned beneath the center of each crate. Several men are required to unload individual components from the crates.
- C. You'll need heavy rubber hammers, good portable drills/screwdrivers, measure tapes (to make sure everything's square) and six to ten 11" locking C-Clamps used to hold beams in place as you insert fasteners. Another tool that will be needed is an 8" – 12" shaft extension for the screwdriver bits (to drive screws in narrow gaps between parts).



Softwall Modular Cleanroom

Components List

The list includes the basic cleanroom components. Quantities vary based on the specifications of each particular unit.

Electrical Service Equipment:


- One (1) master light toggle switch
- One (1) master blower switch (ON/OFF pushbuttons)
- Control boxes (each controls up to 6 lights and 6 fan filter units modules)
- 115 or 230 VAC yellow power cables (normally connected to FFUs and lights before shipping).

PICTURE	USED FOR	DESCRIPTION
	Frame	Black 3/8" X 1/2" Allen Screw Socket Set Cone Point
	Grip Track and Trimming	Screw SS # 12 X 3/4" Phillip Pan Self Drilling
	Support Bar & Joists	Screw SS # 12 X 3/4" Self-Tapping Sheet Metal Screw
	Bit for Trim Strip	Drill Bit , Phillip Power
	Bit for Joist	Drill Bit, 5/16" Magnetic
	Used for Black 3 / 8 Allen Screw	Allen Hex Key
	Ceiling Panels (1-Sided)	Clip 1-Arm Ceiling (If requested)
	Ceiling Panels (2-Sided)	Clip 2-Arm Ceiling (If requested)




Softwall Modular Cleanroom


Cleanroom Frame Components

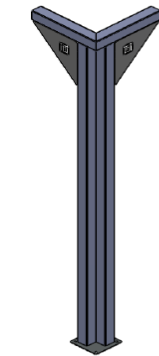


CORNER POST

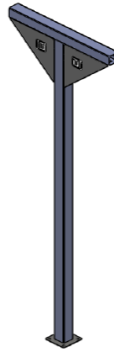

CENTER POST

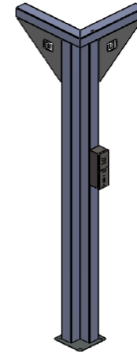


CORNER POWER POST

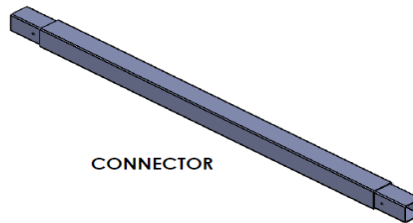


CONNECTOR


Ceiling Components

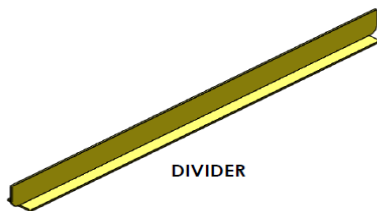


DIVIDER

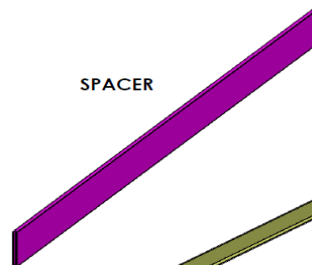

SPACER

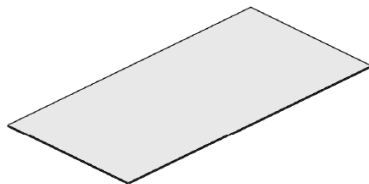

CEILING PANEL

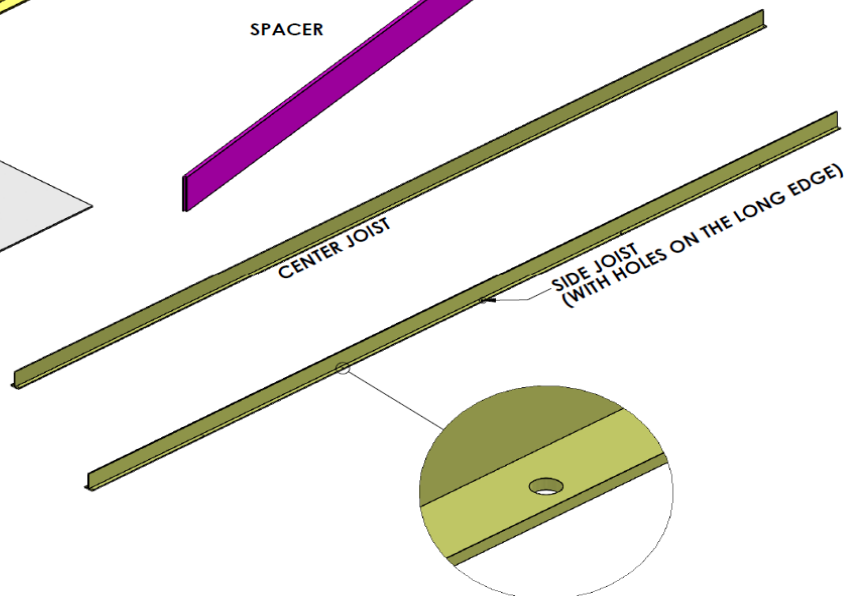

CENTER JOIST


SIDE JOIST
(WITH HOLES ON THE LONG EDGE)




BLACK CLIP



Support Bar Components

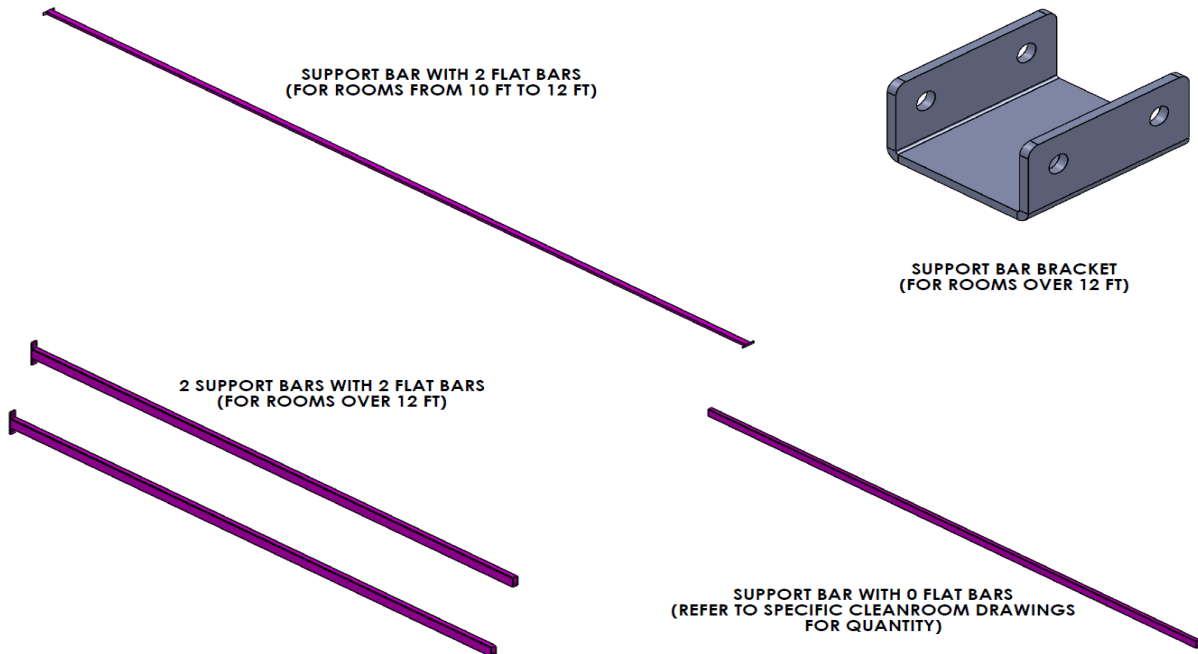


Top Trimming Components

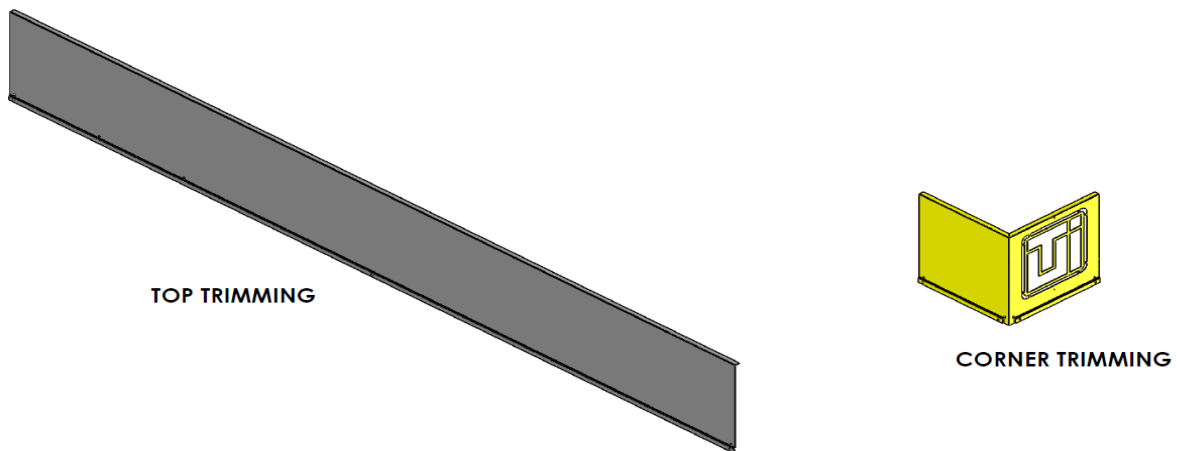



Curtain Panel Components

FLEX-TRACK

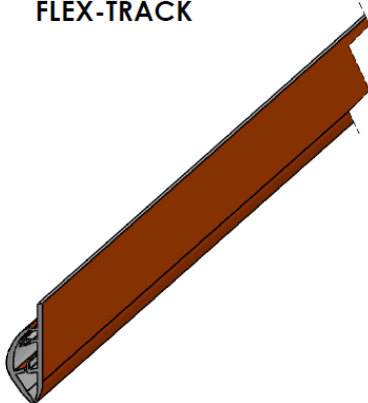

CURTAIN STRIPS



CURTAIN PANELS



Frame Assembly

Installation Note: Before beginning installation of the support frame, familiarize yourself with the placement of the frame members.



IMPORTANT SAFETY NOTICE

Terra Universal Cleanrooms are not designed to support more weight than the blower modules and lighting fixtures originally installed. In particular, the ceiling grid beams are not load bearing and will not support personnel or other additional loads. Placing added weight on the ceiling grid may result in serious damage to the cleanroom and its occupants.


Safety notices supplied by Terra Universal must be affixed at appropriate places on each side of the cleanroom grid.

Prior to shipping, all frame and ceiling members are stamped and then labeled at each end. Refer to the appendix for detailed information on how your cleanroom is numbered.

Before you start assembly, it is mandatory that the floor is level to assure the completed room will fit properly and be rectangular. Failure to level the floor may result in the inability to complete the assembly of the room or the insertion of the blowers, lights or ceiling. For vertical alignment, measure with the leveler before assembly.




Softwall Modular Cleanroom

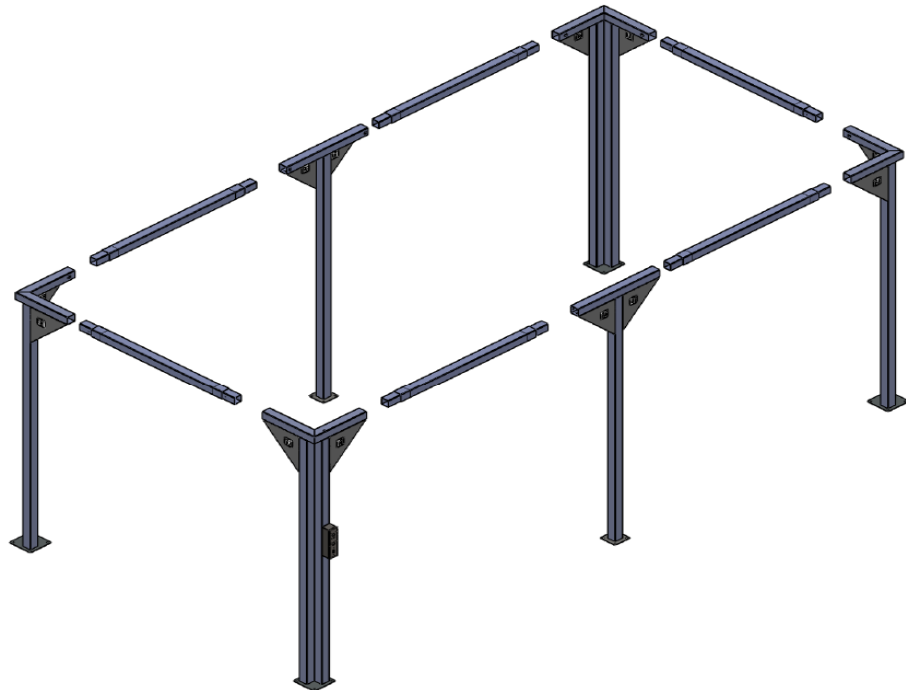

STEP 1: MAKE SURE THE FLOOR IS LEVEL FIRST


STEP 2: REFER TO THE DRAWING TO SEE HOW EACH COMPONENT COMES TOGETHER STARTING AT A CORNER FRAME

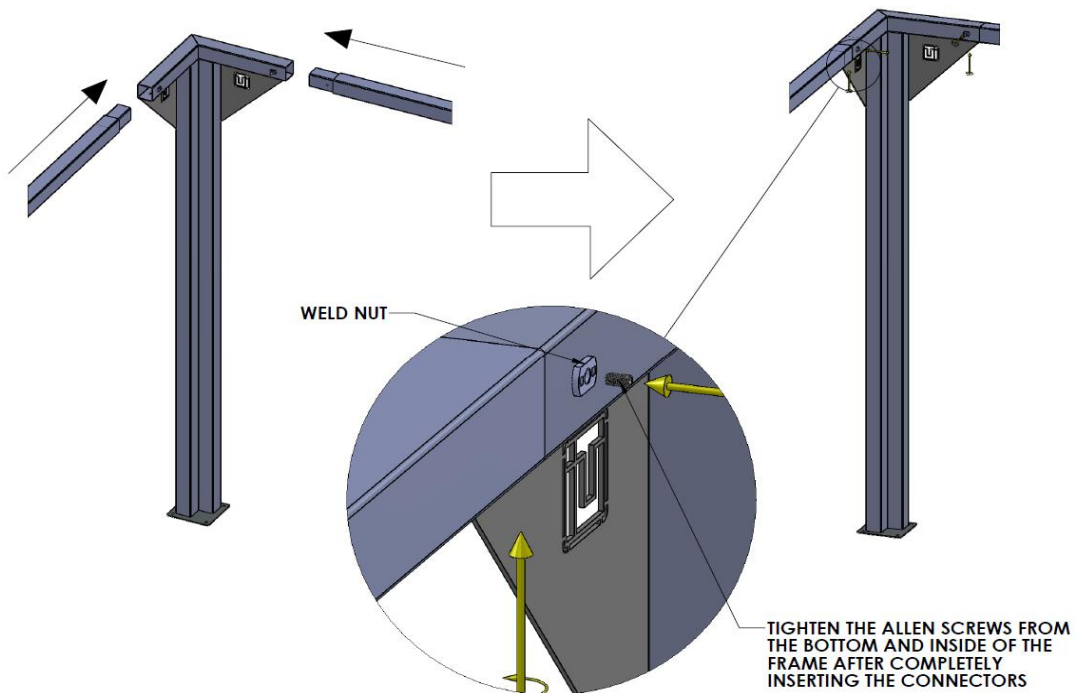

STEP 3



Softwall Modular Cleanroom

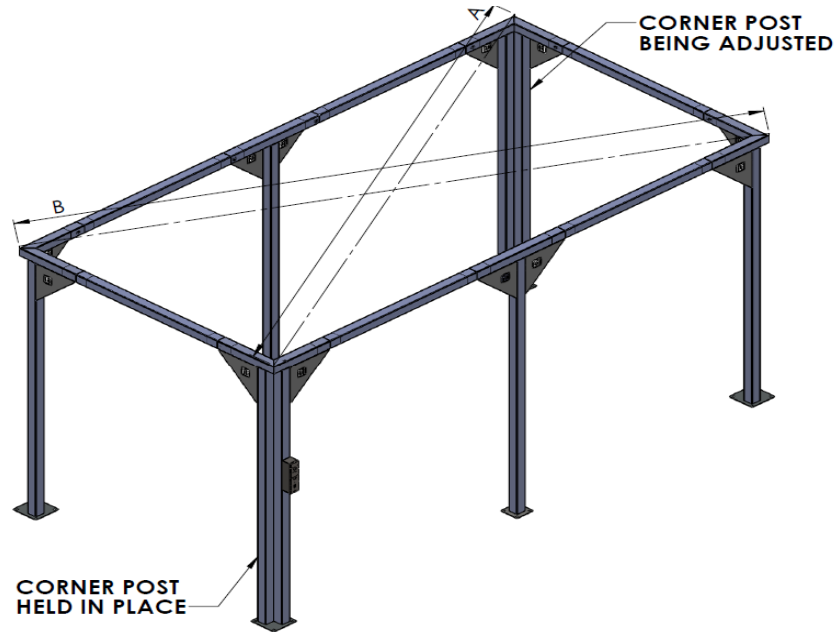

IF THE ROOM IS NOT SQUARED, HAVE SOMEONE HOLD ONE
OF THE CORNERS POSTS DOWN AND ANOTHER PERSON MOVE
THE OPPOSITE CORNER POST EITHER IN OR OUT UNTIL
LENGTH A=LENGTH B

STEP 4

Also, verify that the overall dimensions match the drawing. The frame of the cleanroom can also be anchored to the floor if required.




Softwall Modular Cleanroom

Seismic Anchoring

Because most Softwall Cleanroom models have custom configurations, the number of screws required for seismic anchoring will vary based on the size of the cleanroom. Each post on the frame has a foot with four pre-drilled holes. Be sure to adhere to local seismic codes when anchoring the cleanroom. Terra recommends drilling 6" deep and placing a 3/8" Kwik Bolt TZ (not included) anchor for a minimum embedment of 3-5/8" in each of the foot's holes. This is an example, but local seismic codes may vary.


EXAMPLE DIAGRAM

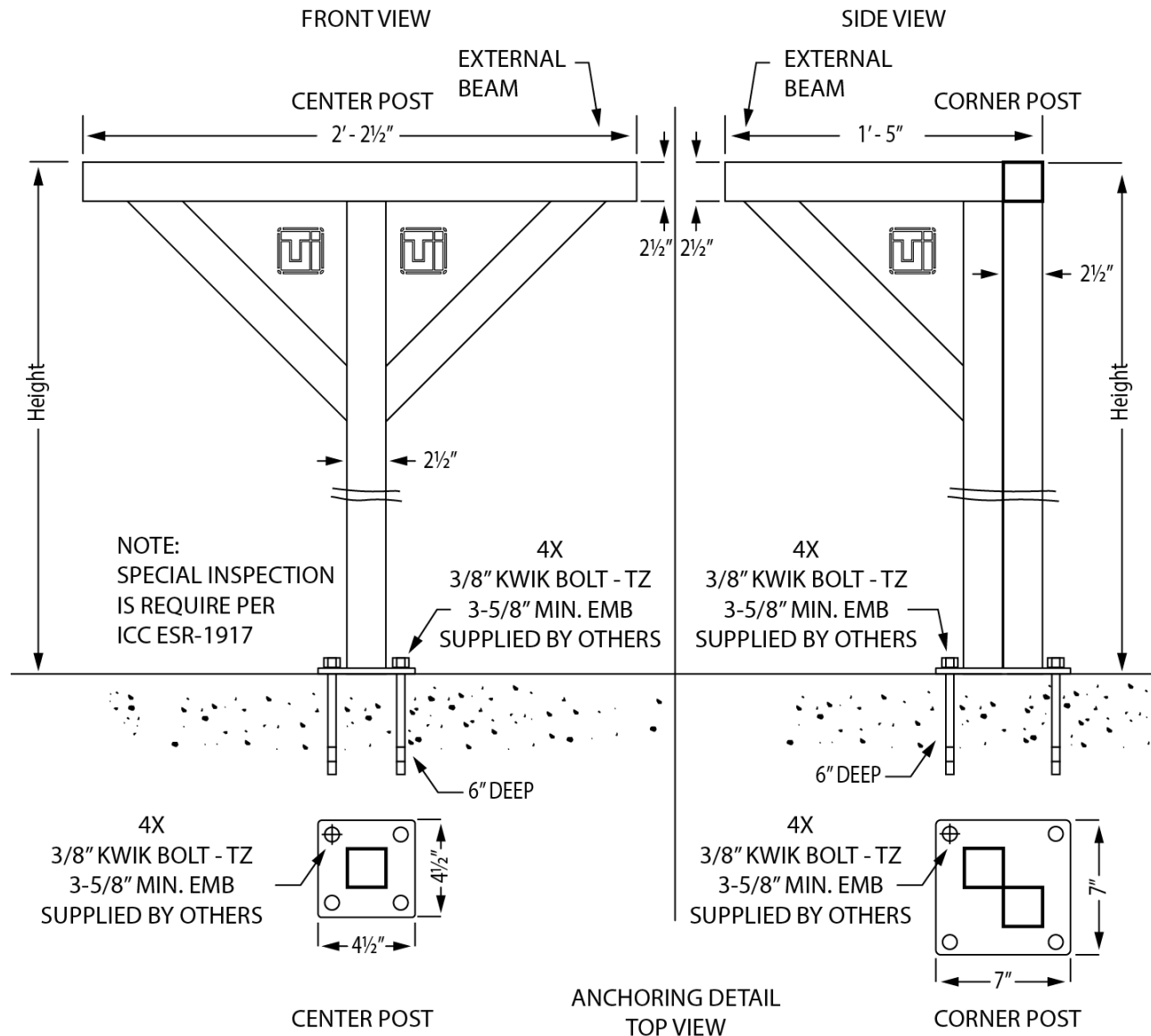



Softwall Modular Cleanroom


Joist and Divider Installation

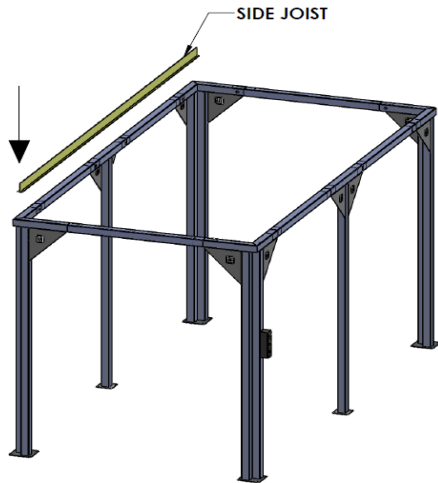

JOISTS ARE BOWED TO COMPENSATE FOR THE WEIGHT OF ALL THE FFUS, LIGHTS, PDMS, AND OTHER CEILING ACCESSORIES


NOTE: ONLY SIDE JOISTS HAVE EXTRA FASTENING HOLES

STEP 1

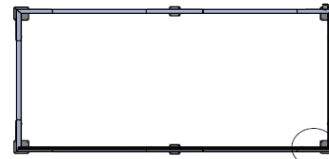


TOP VIEW

THERE ARE NO PREDRILLED HOLES ON THE FRAME

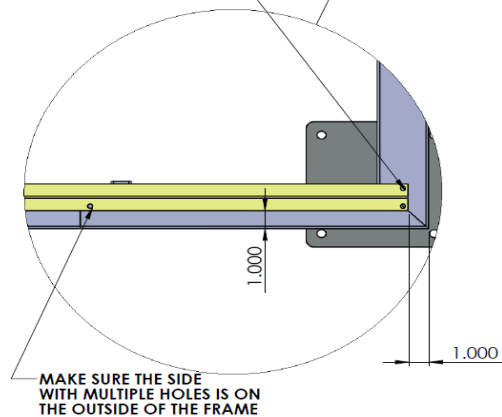


STEP 2

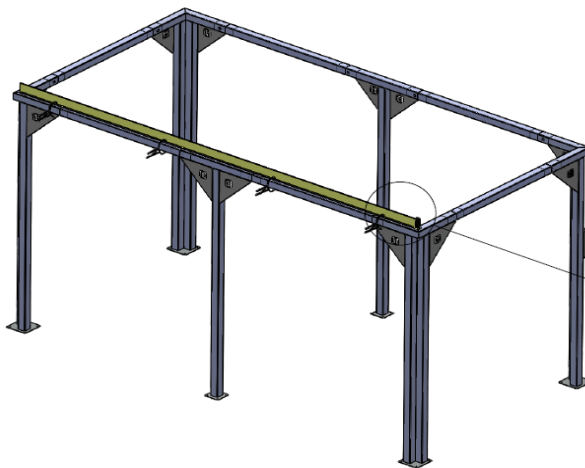

USE SELF TAPPING SHEET METAL SCREWS TO DRILL A HOLE AND ATTACH THE SIDE JOIST TO THE FRAME

CLAMP THE JOIST TO THE FRAME TO SCREW IT ON


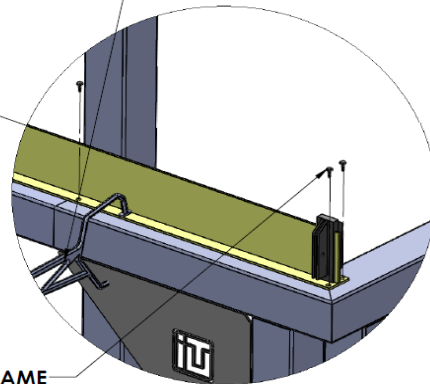


REMOVE THE CLAMPS AFTER SCREWING THE JOIST TO THE FRAME



STEP 3



Softwall Modular Cleanroom

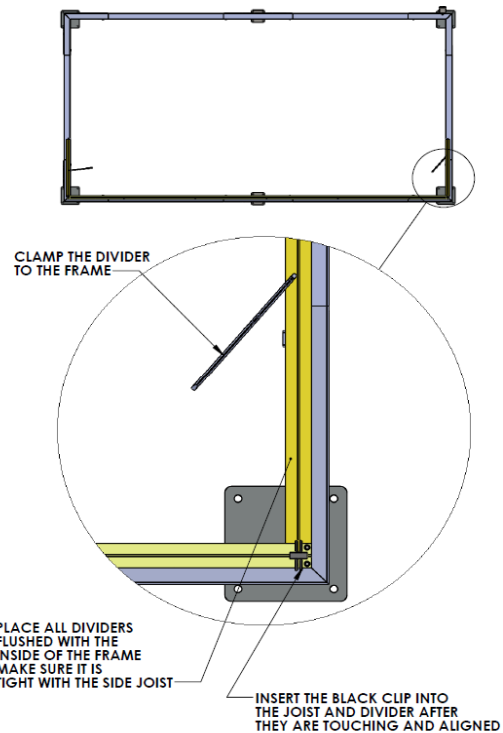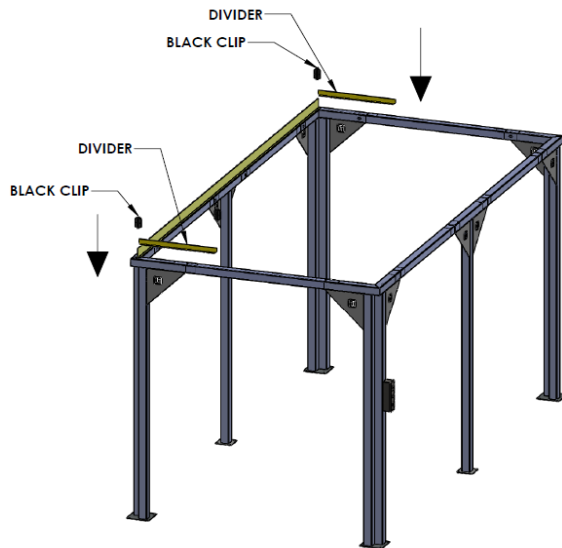

STEP 4

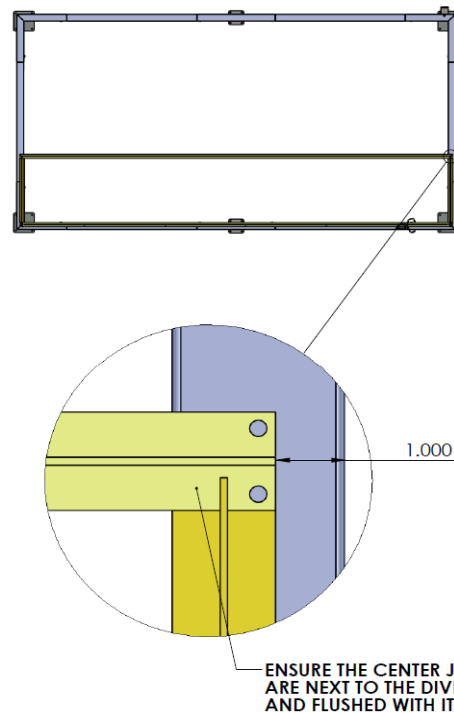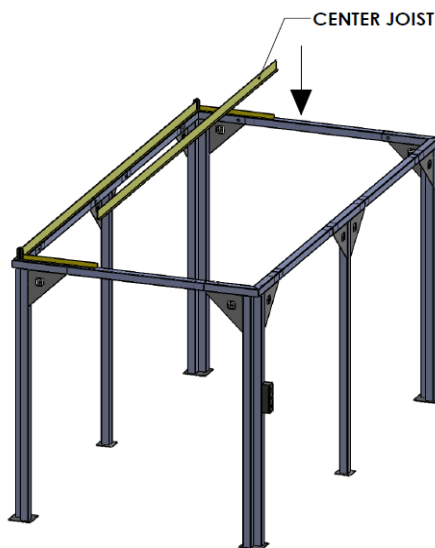


STEP 5


NOTE: CENTER JOISTS DO NOT HAVE HOLES ALONG THE LONG EDGE



Softwall Modular Cleanroom

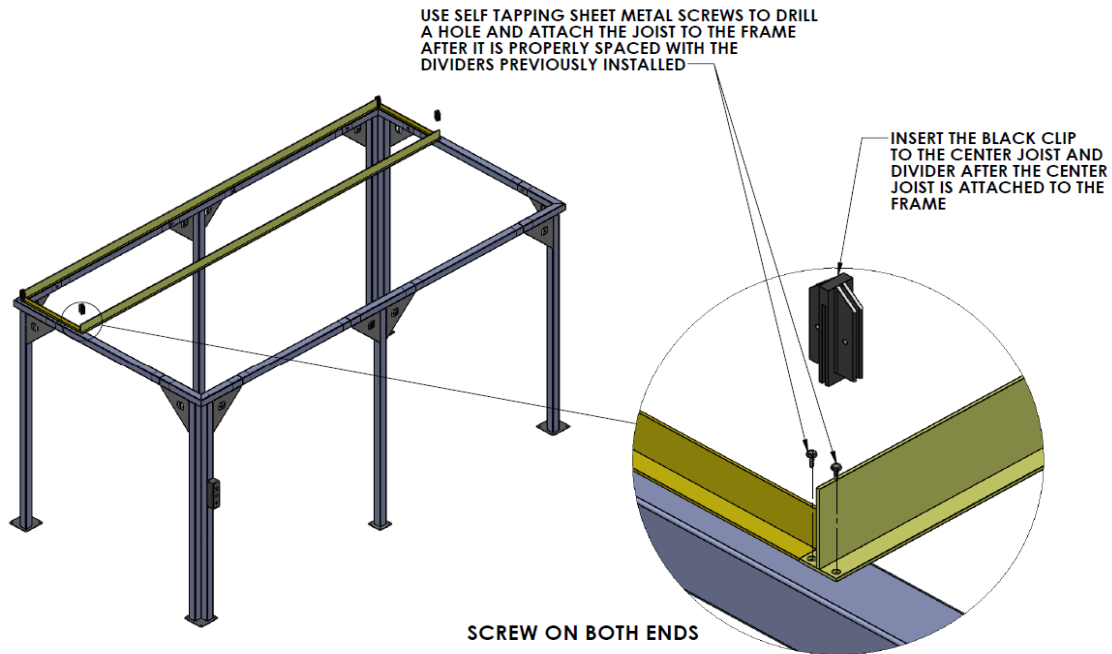



STEP 6

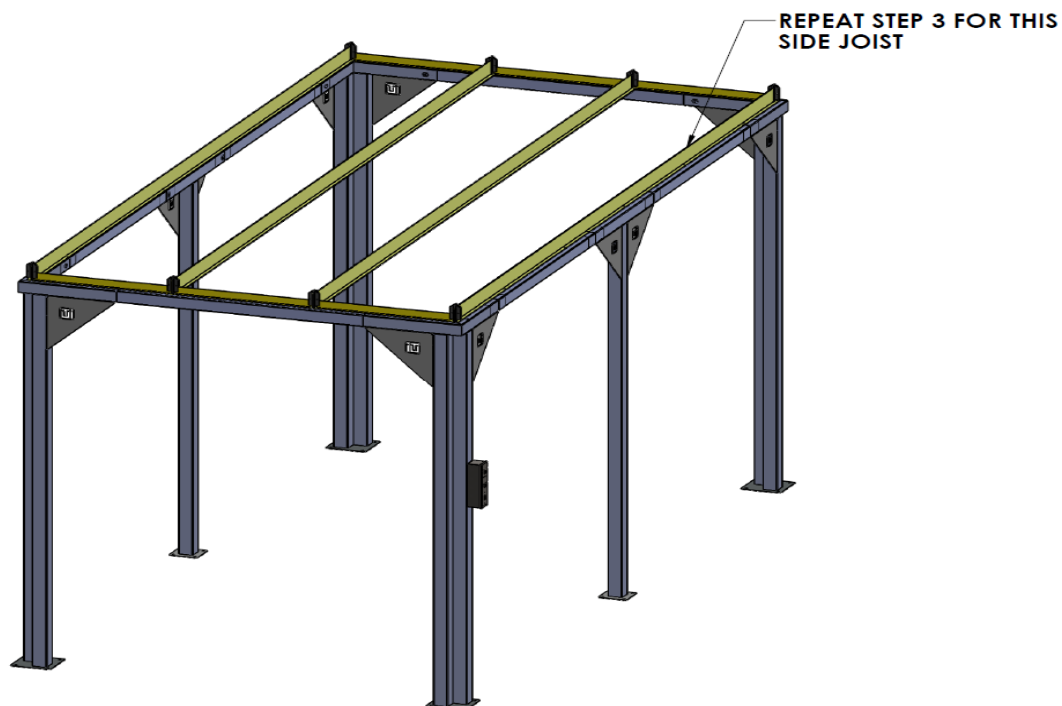



Softwall Modular Cleanroom

Use of the Spacer

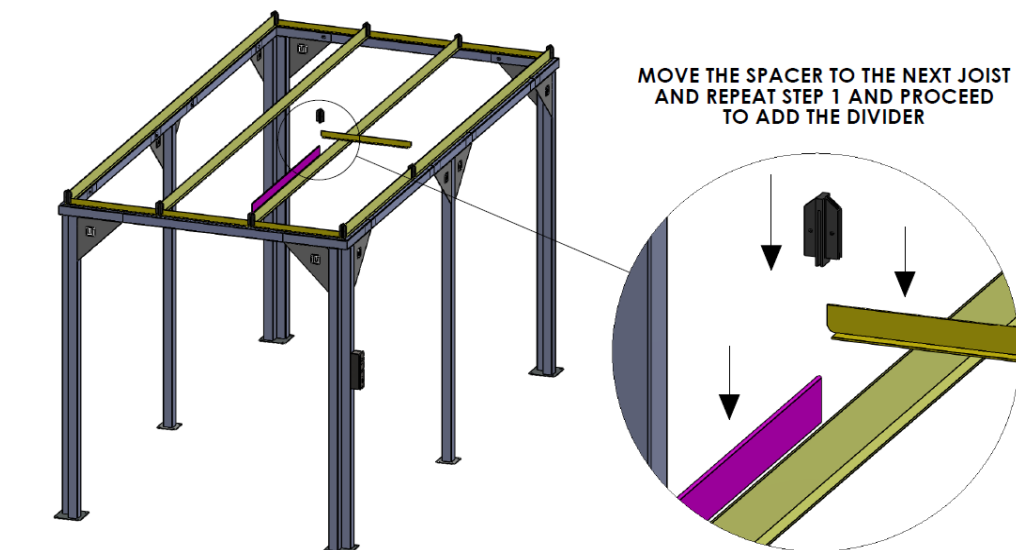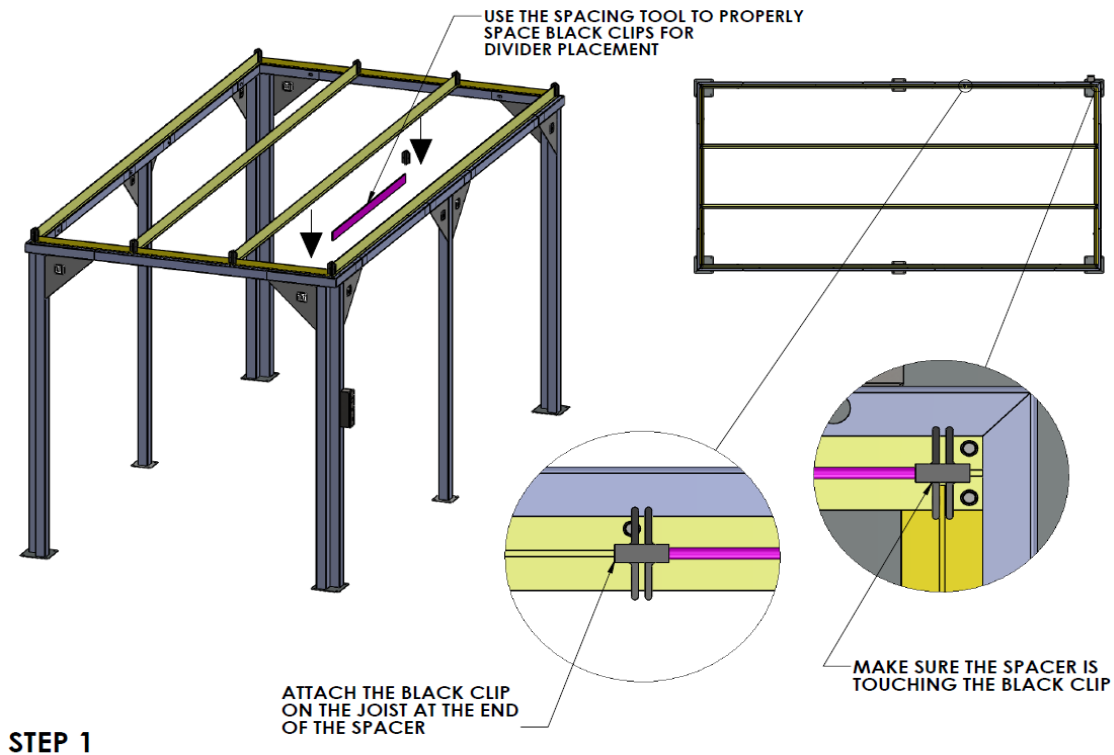



Softwall Modular Cleanroom



STEP 3


**REPEAT STEPS 1-2 UNTIL THE
CEILING GRID IS COMPLETE**

Recheck the frame to ensure that the structure is square.




Softwall Modular Cleanroom

Support Bar Installation (For Cleanrooms from 10 ft. to 12 ft.)

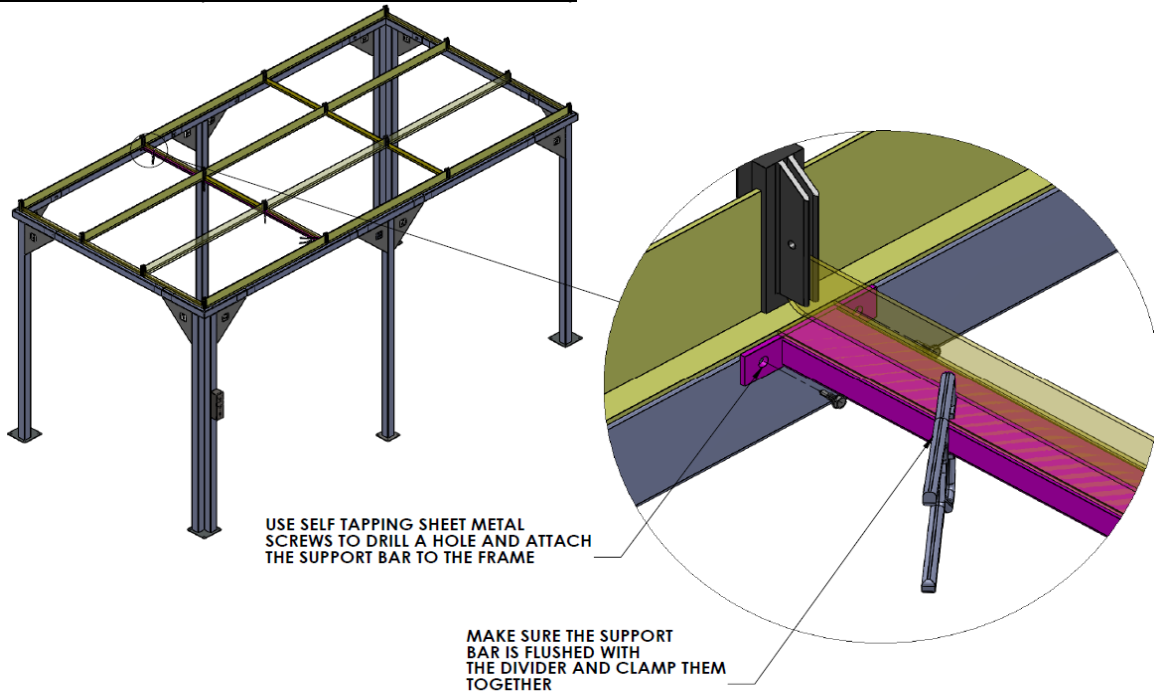

STEP 1


REPEAT THIS STEP FOR THE OPPOSITE END OF THE ROOM

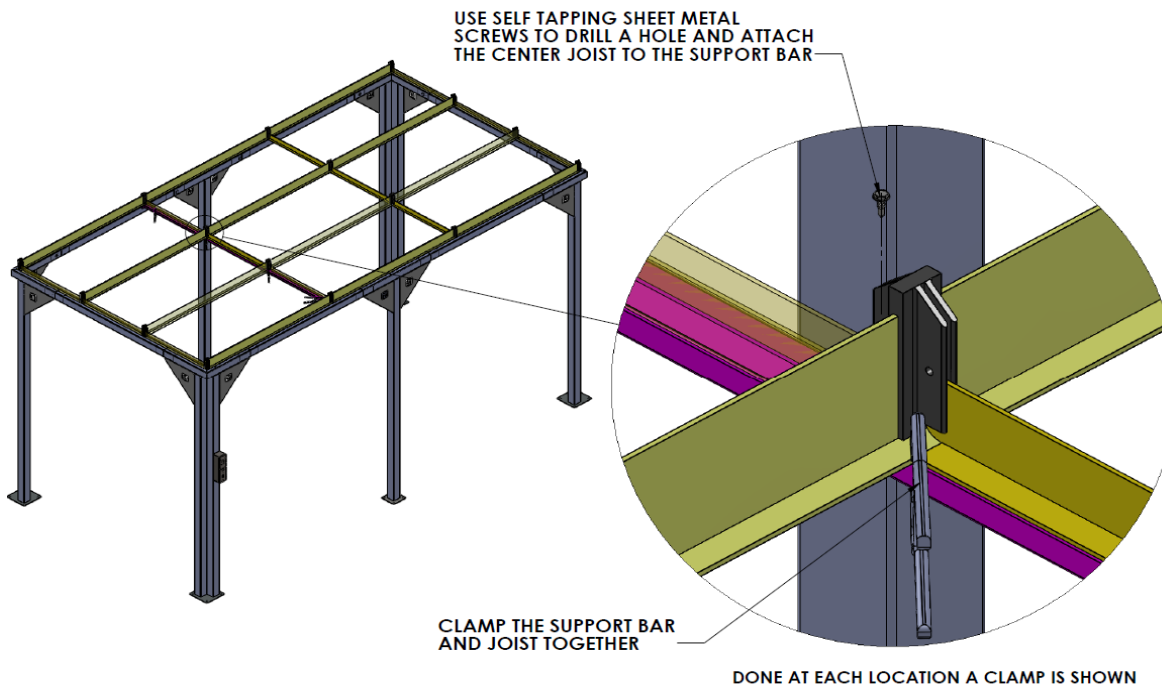


STEP 2




Softwall Modular Cleanroom

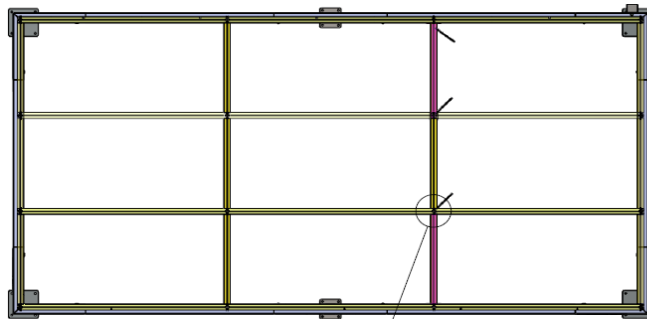

TOP VIEW

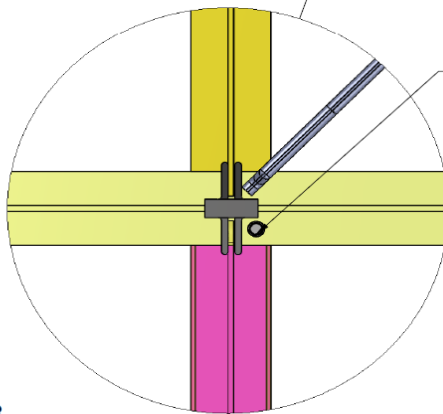

USE SELF TAPPING SHEET METAL
SCREWS TO DRILL A HOLE AND
ATTACH SUPPORT BAR TO THE JOIST

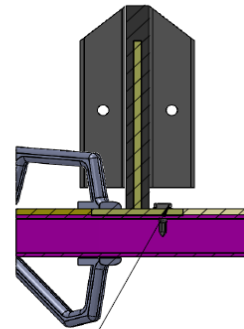


MAKE SURE THE SCREW GOES
THROUGH THE JOIST AND SCREWS
ONTO THE SUPPORT BAR.

STEP 3




Softwall Modular Cleanroom

Installing 2 Support Bars (For Rooms Bigger Than 12 ft.)

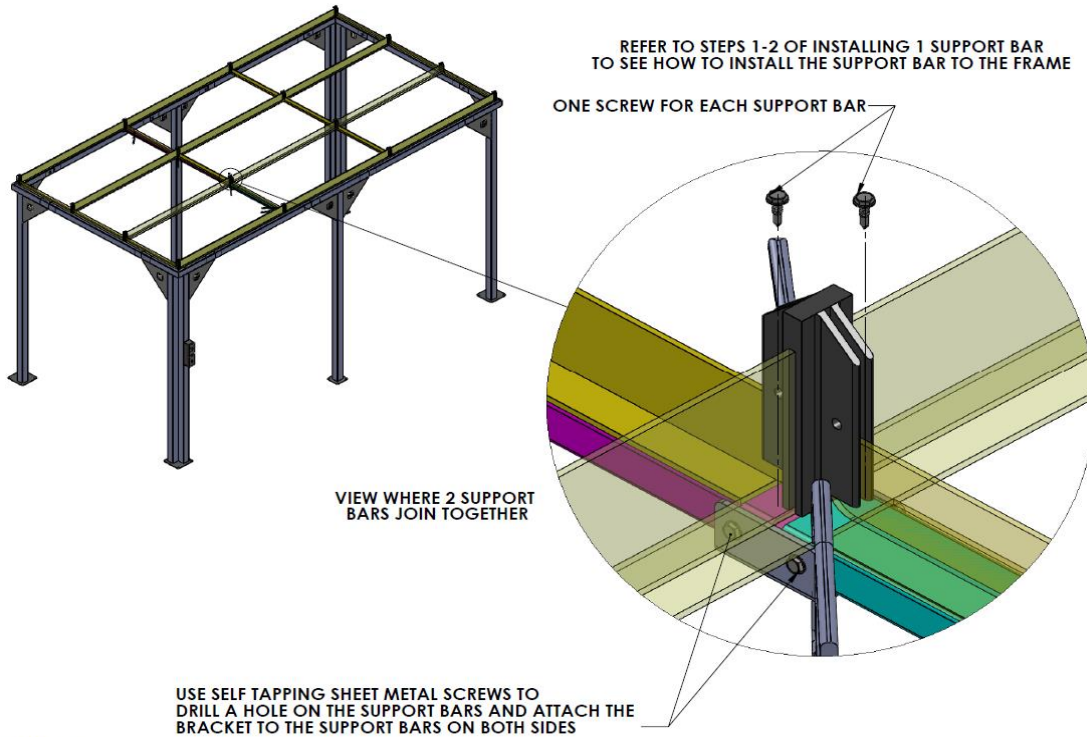

STEP 1

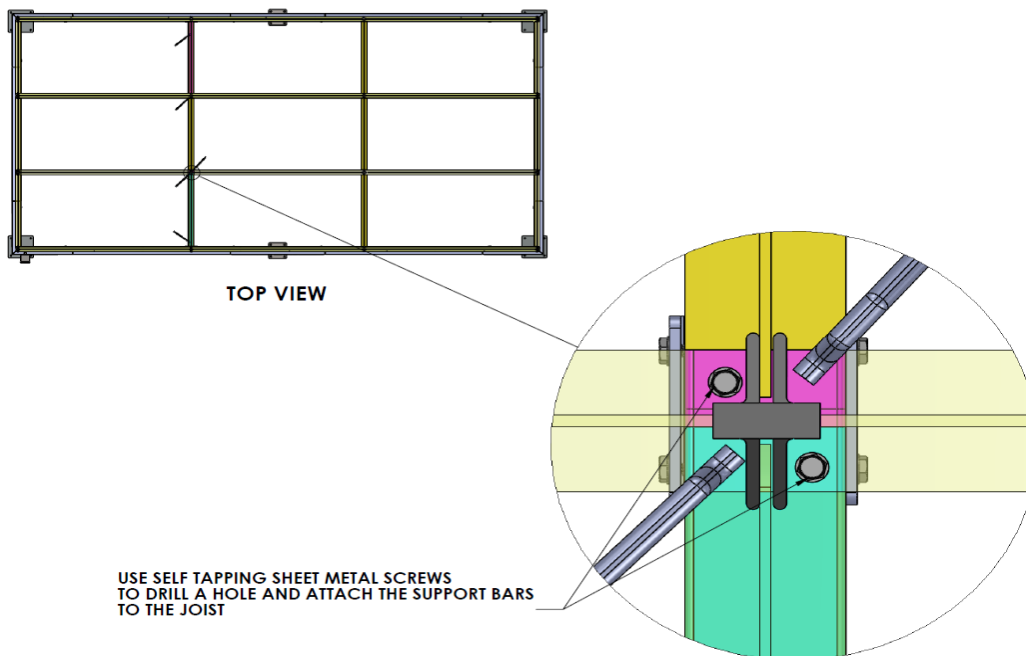

STEP 2


PAGE 26 OF



Softwall Modular Cleanroom


© Copyright 2021 Terra Universal Inc. All rights reserved. • Revised

Top Trimming Installation

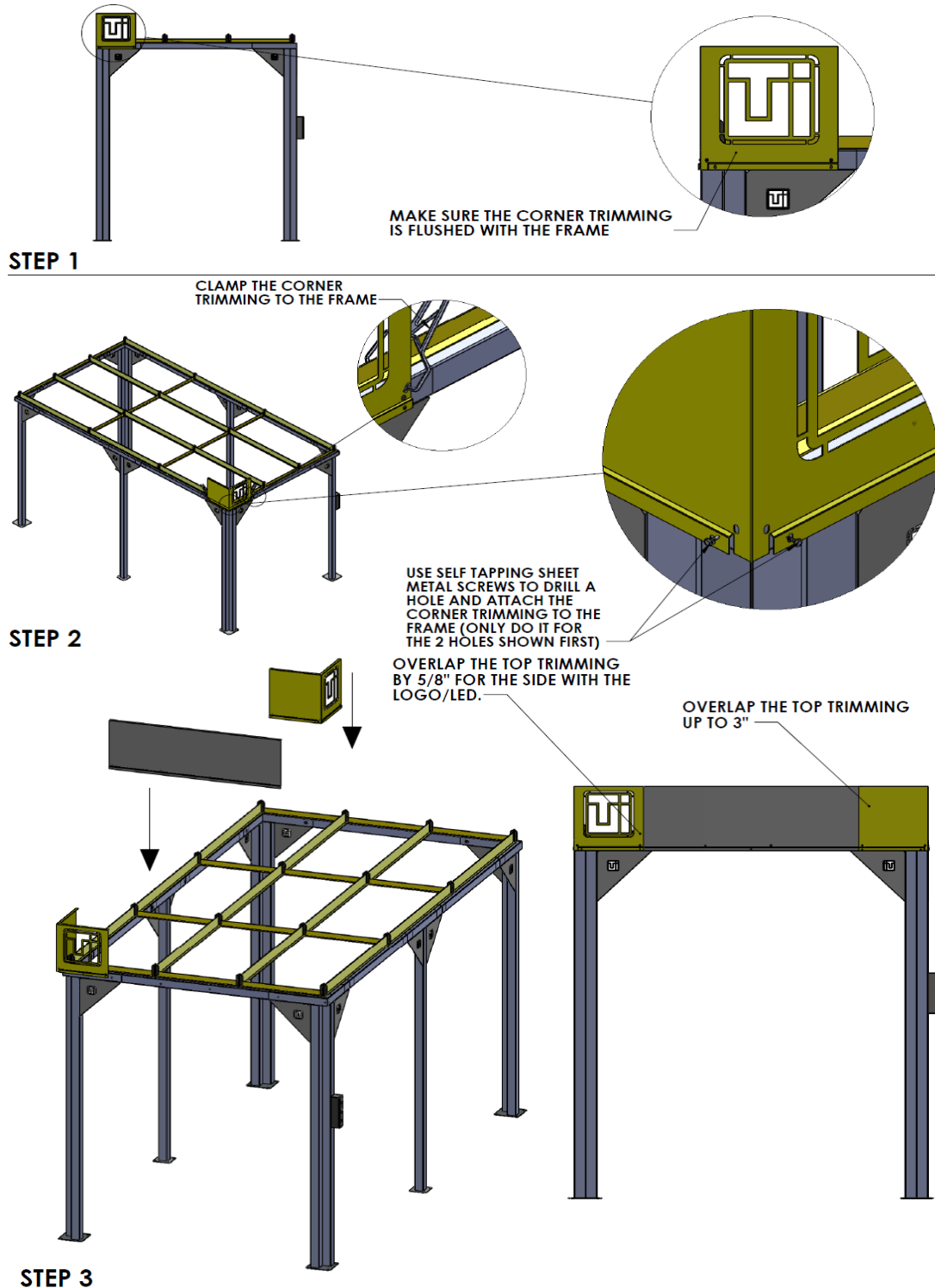




Softwall Modular Cleanroom

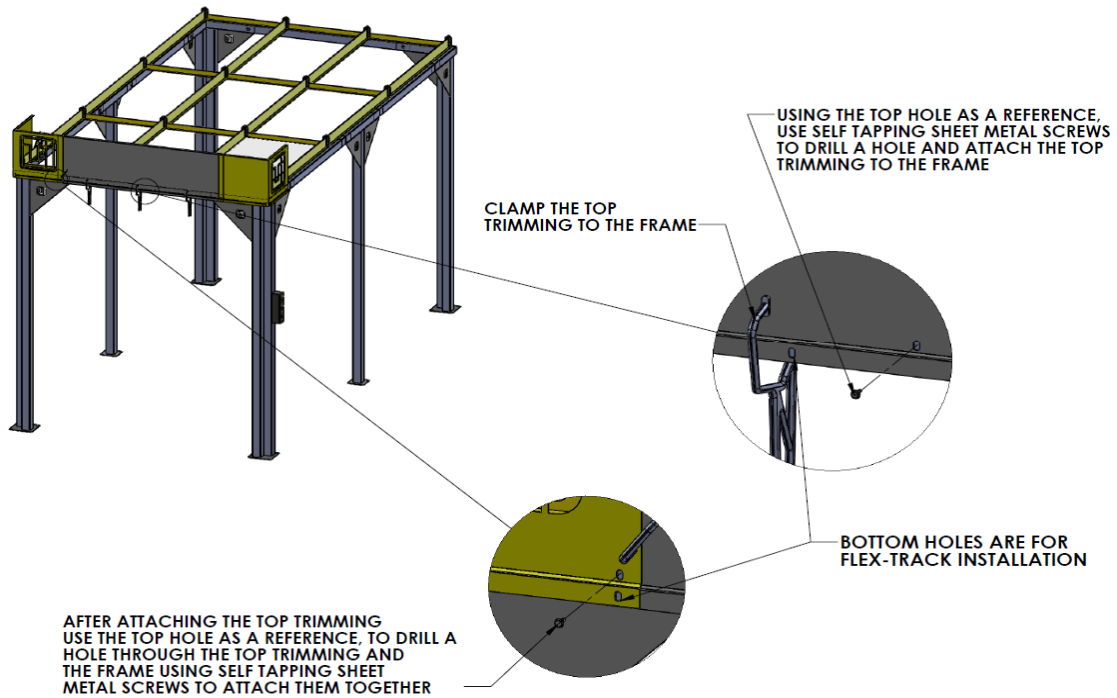


STEP 4

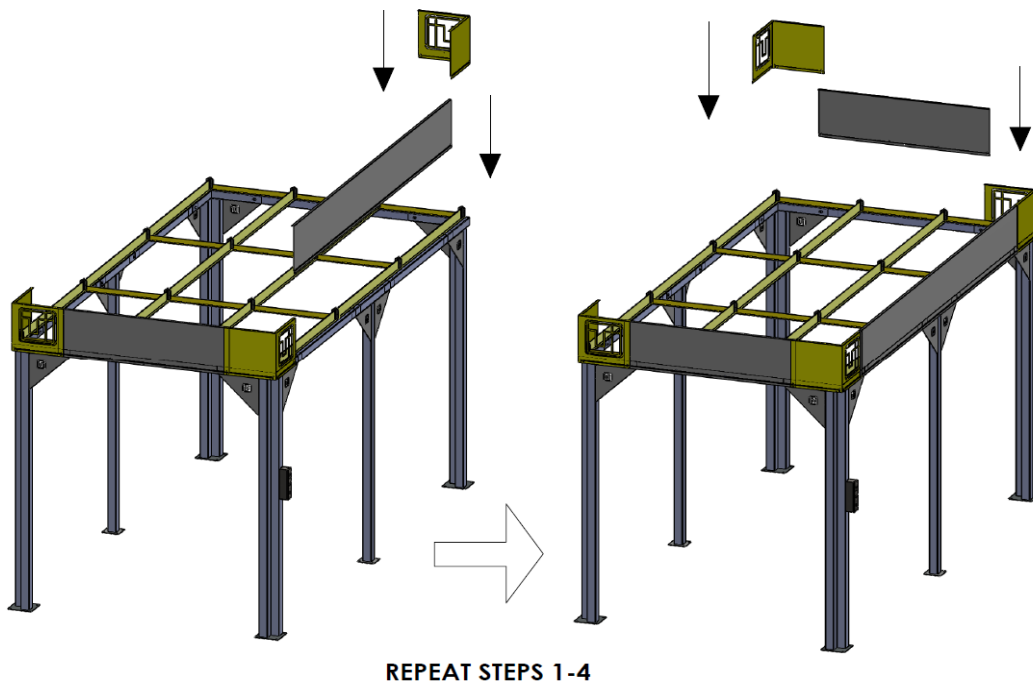

STEP 5



Softwall Modular Cleanroom

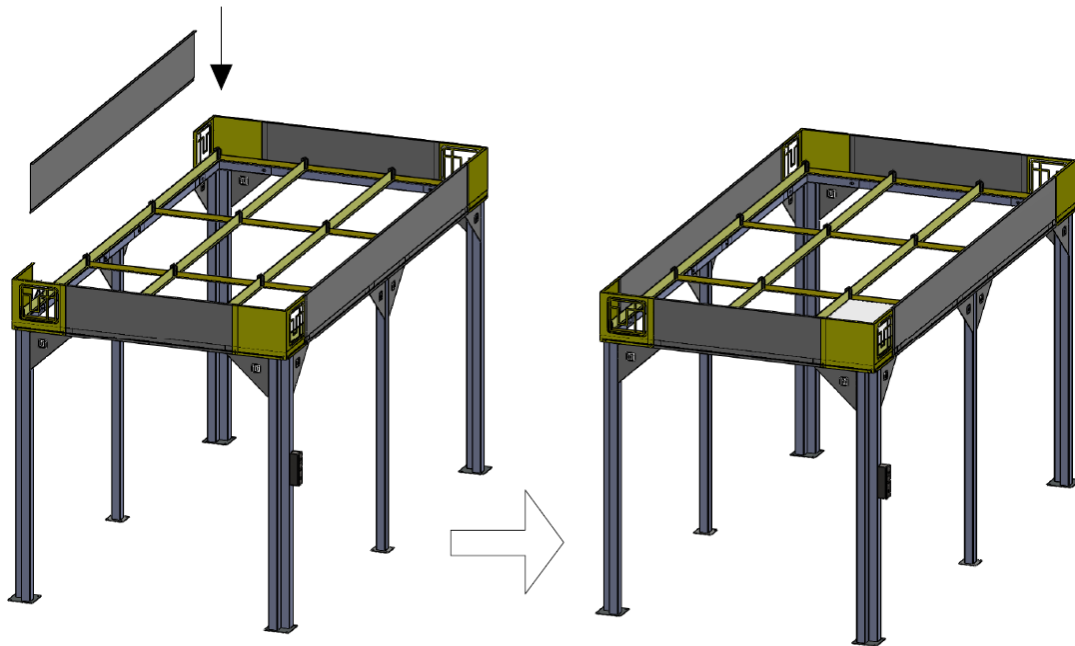


REPEAT STEPS 3-4

STEP 6

Note: To prevent scratching, leave protective plastic on steel trim until assembly is complete.

Installing the Ceiling Modules

**DROP CEILING ACCESSORIES
ONTO THE OPENINGS**

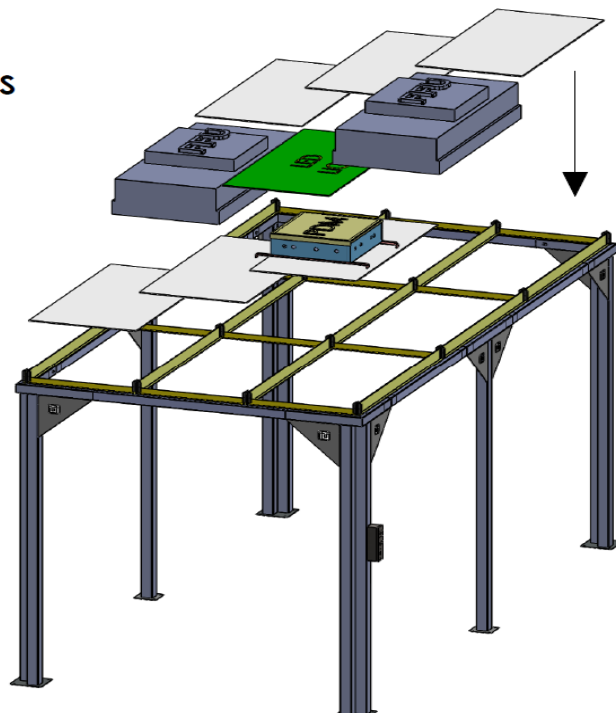

**NOTE: REFER TO THE ELECTRICAL
SECTION OF THE MANUAL FOR
WIRING DETAILS**



Softwall Modular Cleanroom

If provided, secure down with Terra-provided Cleanroom Ceiling Clips. Use the 1-Arm Clip for joists with only one adjacent ceiling panel and use the 2-Arm Clip for joists with two adjacent ceiling panels. See Figures 1 thru 4.

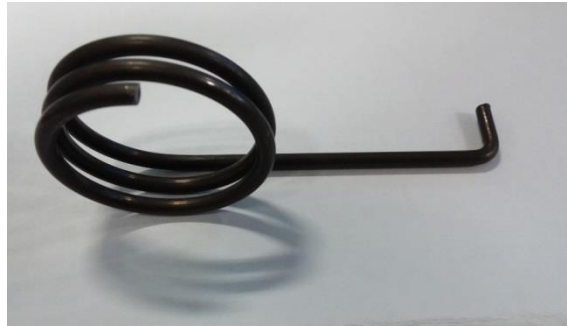


Figure 1: 1-Arm Ceiling Clip



Figure 2: 2-Arm Ceiling Clip



Figure 3: 1-Arm Ceiling Clip holds down a ceiling panel with a FFU installed on the opposite side of the joist.


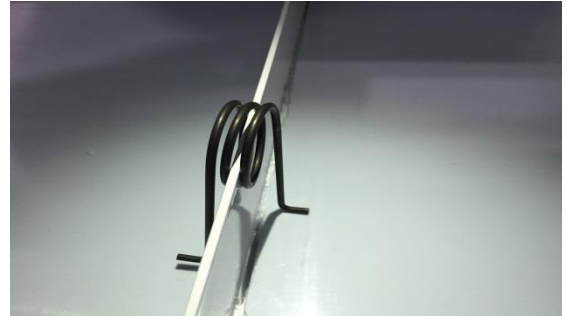


Figure 4: 2-Arm Ceiling Clip simultaneously holds down two ceiling panels.




Softwall Modular Cleanroom

Curtain Panel Installation

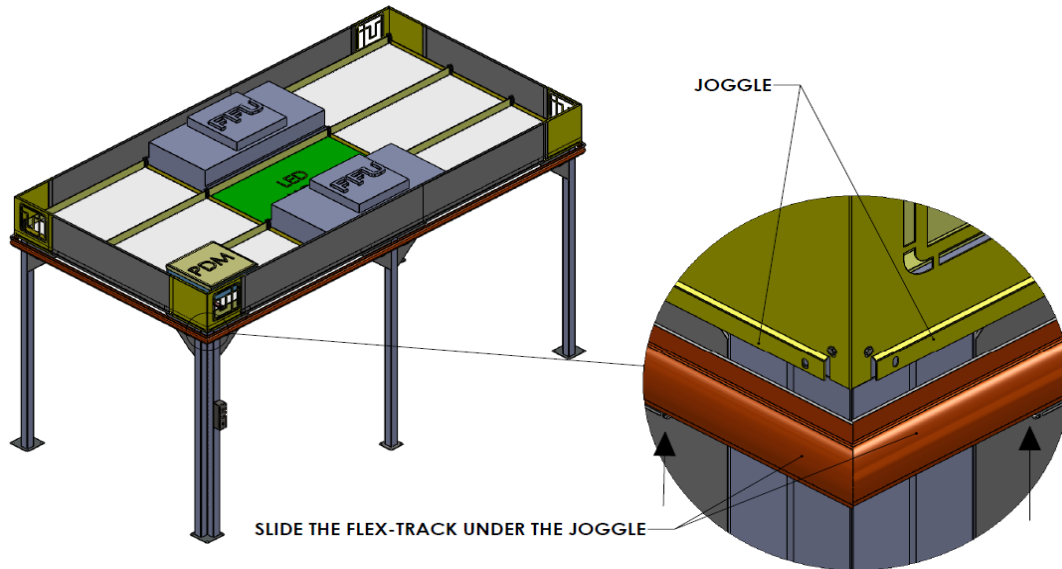

MAKE SURE THE FLEX-TRACKS TOUCH AT THE CORNERS AND TRIM THE FLEX-TRACK IF NECESSARY


STEP 1

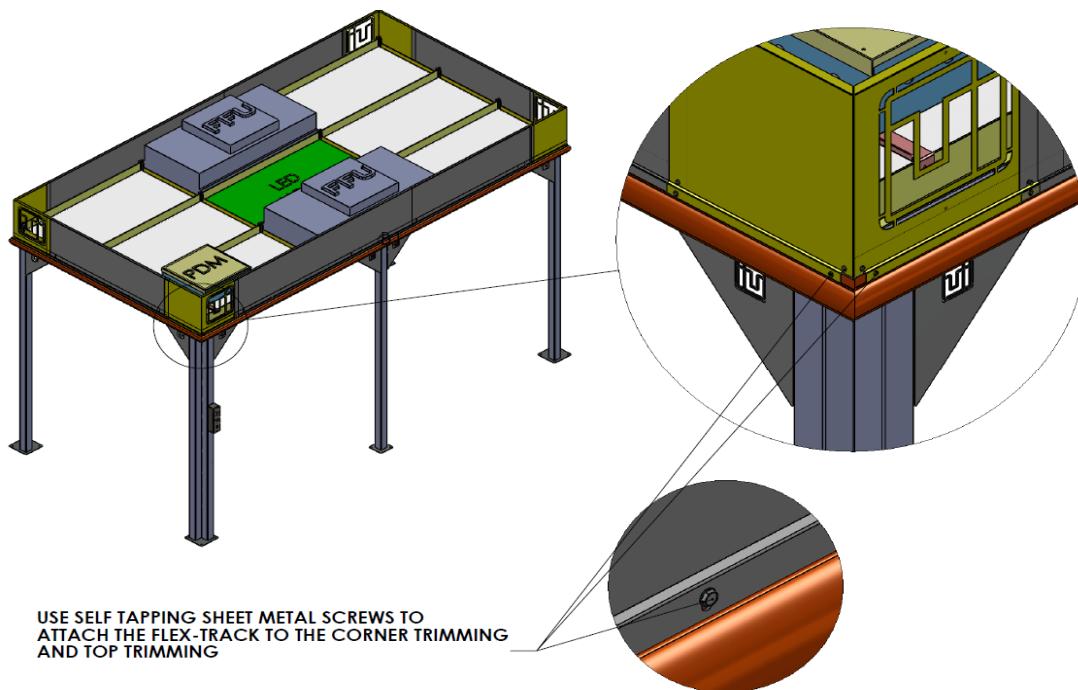

STEP 2




Softwall Modular Cleanroom

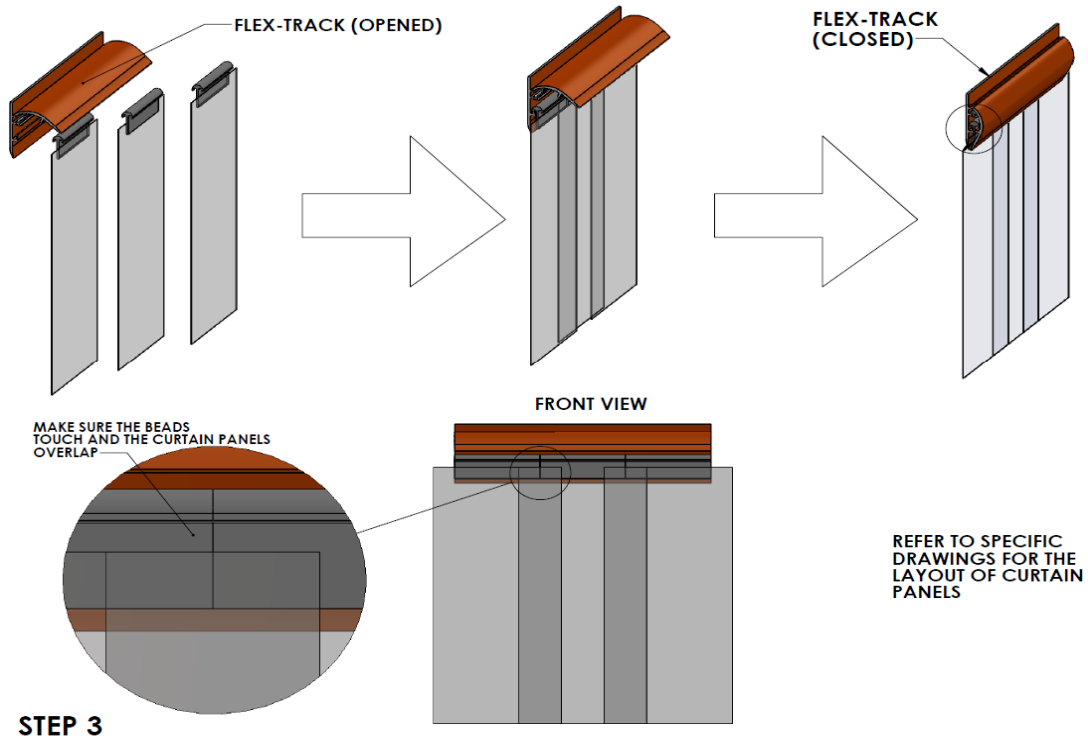



3.0 Electrical Wiring



Installing the PDM

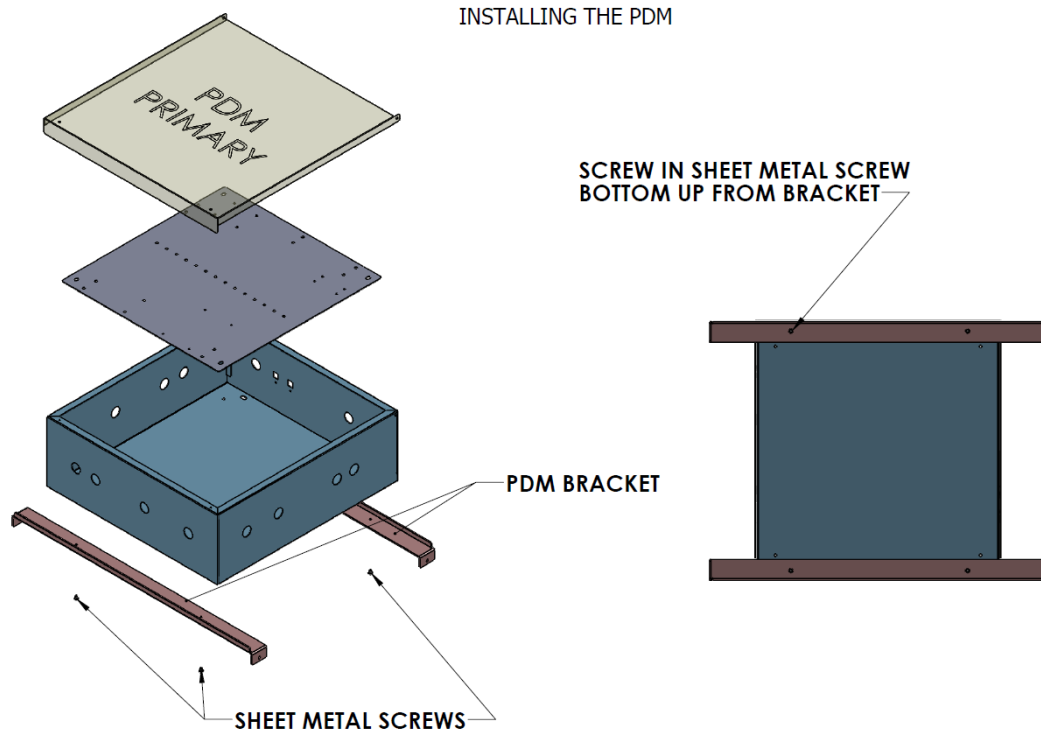




Softwall Modular Cleanroom

REFER TO THE DRAWING
TO DETERMINE WHERE TO
PUT THE PDM

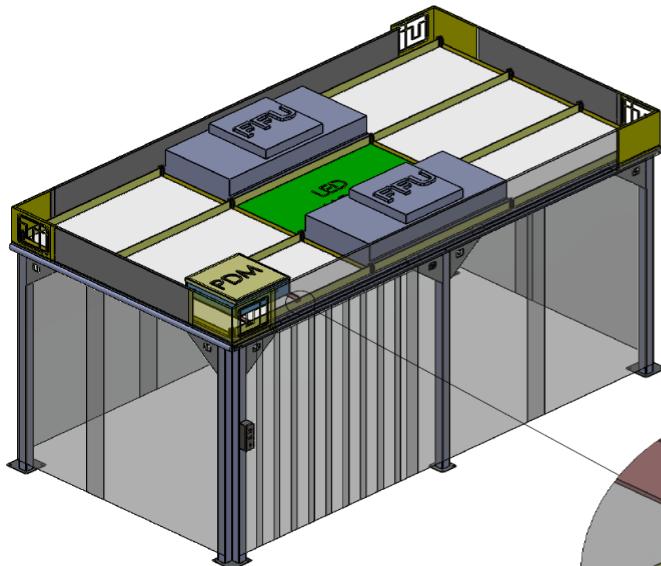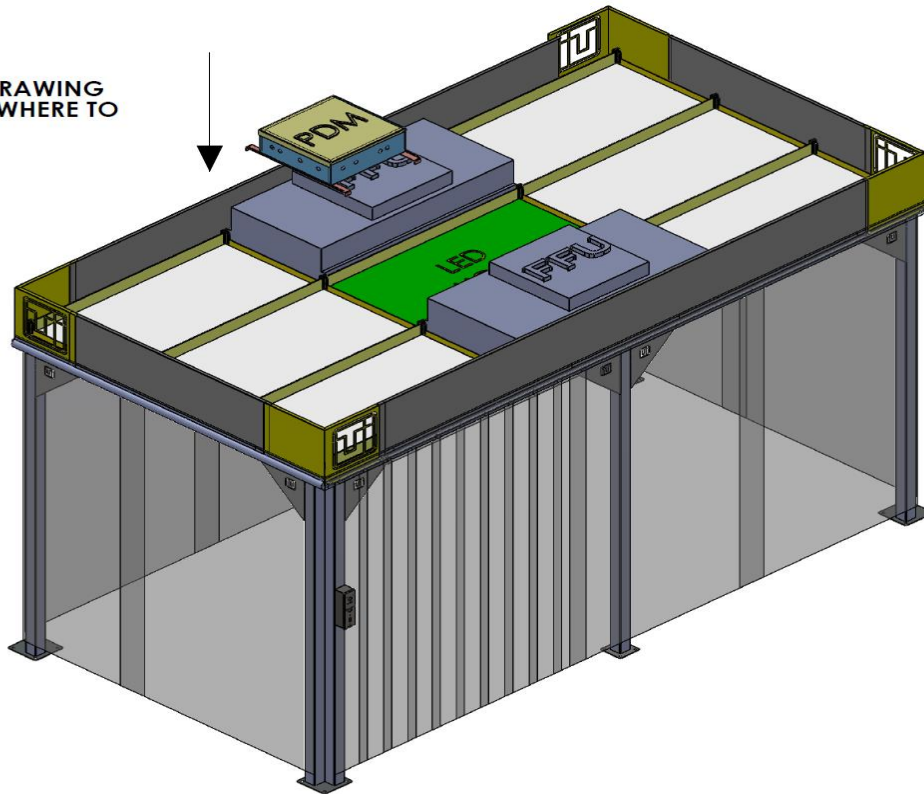

USE SELF TAPPING SHEET METAL SCREW
TO DRILL A HOLE ON THE JOIST TO
ATTACH THE PDM

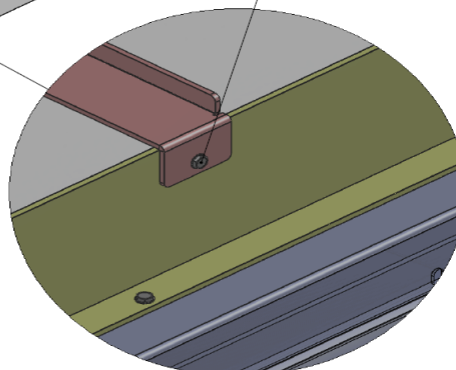



Softwall Modular Cleanroom

Use the provided yellow power cables to connect the components according to the labels on the PDM. 4-pin yellow cables are for the FFUs and 3-pin yellow cables are for the light fixtures.



Figure 1: PDM connection for the Control Panel



Figure 2: PDM connection for light fixtures



Figure 3: PDM connections for the fan/filter units



Figure 4: Cord grip on fan/filter unit housing



Operation Status Indicator Lights Wiring

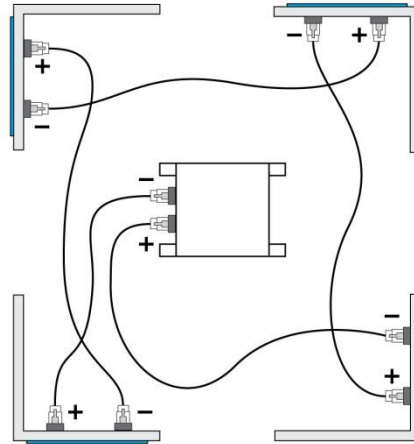


Figure 5: Wiring Schematic for Operation Status Indicator Lights.

Use the provided low-voltage phone cable for connections

To test for proper operation, hard wire the unit to a grounded 115VAC/60Hz (or 220VAC/50Hz, where applicable) power source and turn the switches ON. All electrical modules should come on.



Fan Filter Unit Hard Wiring

! WARNING: Disconnect the unit from the electrical power source before attempting any service.

To preserve the UL listing of Terra Fan Filter Units, Terra must configure them for hard-wiring. Cords are dressed to simplify this operation, which typically does not require an electrician.

1. Remove the blue prefilter, which rests on top of the fan filter unit (Figure 6).
2. Unscrew metal top panel of the electrical housing (Figure 7) to expose wiring junction (Figure 8).
3. Install cord grip as shown (Figures 9, 10) into appropriate opening.
4. Thread the AC power chord through the cord grip and connect to wire port as shown (Figure 11). Match the colors of the wiring, as indicated on the charts below. Tighten retaining screws to fasten wires.
5. Tighten cord grip to secure AC power cord.
6. Replace the metal panel on top of the electrical housing.
7. Replace the blue prefilter.



Fig 6



Fig 7



Fig 8



Fig 9



Fig 10



Fig 11

120V FFU Wiring				
Cable Wire:	White	Black	Red	Green
Function:	Neutral	Line	None (cut)	Ground
FFU Connector:	White	Black	None	Green

220V FFU Wiring				
Cable Wire:	White	Black	Red	Green
Function:	None (cut)	Neutral	Line	Ground
FFU Connector:	None	Black	Red	Green




Softwall Modular Cleanroom

Fan Filter Unit Speed Configuration with PDM





To avoid thermal overload on FFU's with Permanent Split Capacitor (PSC) motors, do not use LOW speed switch when the FFU is connected to a PDM.

The electric circuit design of the PDM limits speed configurations of FFU's with PSC motors. All FFU's with PSC motors should be switched to MEDIUM or HIGH speed setting when connected to a PDM. Thermal overload can occur when one FFU is set to LOW, and the others are set to MEDIUM or HIGH speed settings. The FFU motor set to LOW can draw excess current from the PDM which can lead to thermal overload on the motor, motor failure, and/or electrical hazards.

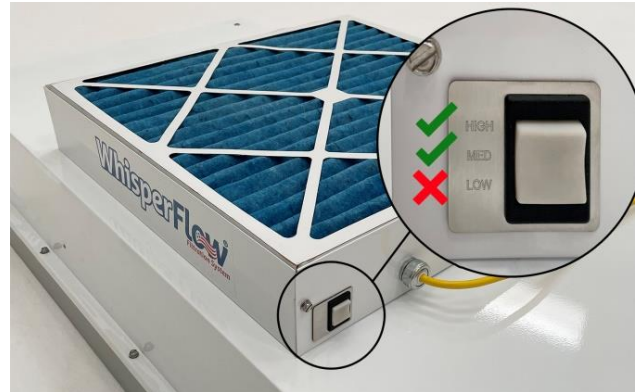


NOTES:

1. Setting all FFU's with PSC motors to LOW does not cause an immediate problem. However, it is not advised as there is a risk of accidentally adjusting one FFU's speed setting to HIGH or MEDIUM, which could lead to thermal overload on a FFU motor that is still on LOW speed.
2. "Night service" mode from the control panel is unaffected and operates correctly when all individual FFU's are switched to HIGH or MEDIUM speed settings.
3. This limitation does not apply if the FFU with PSC motor is used individually without a PDM, and powered directly from a standard wall outlet.
4. This limitation does not apply to Smart® FFU's with Electronically Commutated (EC) motors. Smart® FFU's do not use PSC motors.

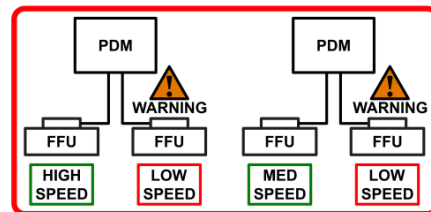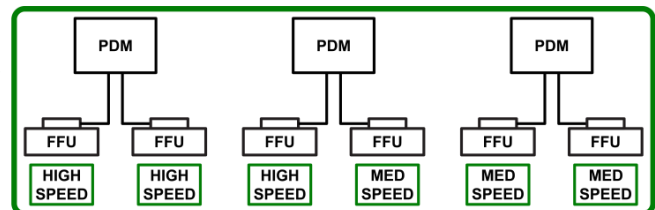

SEE NOTE #1

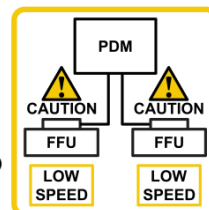



4.0 Operation and Maintenance

Operation of the Modular Cleanroom is controlled by an on/off switch for lights and push buttons for the fan filter units, located near the access door. The Modular Cleanroom typically requires minimal maintenance. The filters provide effective operation for years under typical operating conditions. In fact, filter efficiency increases as the filter captures more and more particles. The filter does not require replacement until the backpressure it generates increases to the point that the system can no longer provide an adequate airflow velocity to maintain required particle counts. To monitor this condition, periodic testing with a particle counter and air velocity meter is recommended.


Filter Replacement

To replace the filter, disconnect system power and remove the fan filter unit from the top of the structure. Then position it to come through the 2' x 4' opening and into the cleanroom. See chart below for ordering information. Detailed replacement instructions on the next page.



WARNING: Never attempt to service or replace the filter without first disconnecting system power to the fan filter unit.

Terra Universal FFU Filter		
Type	Filter Size	TUI Part #
HEPA	2 X 2	6601-27
	2 X 3	6601-26
	2 X 4	6601-25
ULPA	2 X 2	6601-30
	2 X 3	6601-29
	2 X 4	6601-28

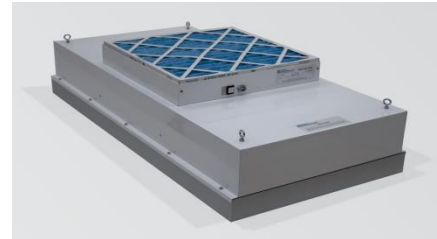

Removal and Replacement of the HEPA/ULPA Filter (Standard Unit)



Warning: Disconnect the unit from the electrical power source before attempting any service.



Warning: The Standard Filter is protected with an expanded metal face screen. This is never to be used to handle the filter. It is only for protection against an accidental touch of the filter. Handle filter only by the frame.



Softwall Modular Cleanroom

Tools Required: Phillips Head Driver

Step 1: Disconnect yellow power cable and remove unit from ceiling (see next page for recommended procedure).

Step 2: Remove the 10 screws holding the HEPA/ULPA filter to the lid assembly.


Step 3: Lift the lid assembly off the HEPA/ULPA filter (see figure below). Discard the used filter as per applicable regulations.

Step 4: Carefully attach the new filter, being sure not to touch or otherwise damage the filter face.

Step 5: Lift out the old pre-filter and drop in the new one.

Step 6: Position the unit back in the ceiling grid and reconnect the unit to its power supply.

Note: Before replacing with a new HEPA/ULPA filter, carefully inspect the new filter for any visible damage.

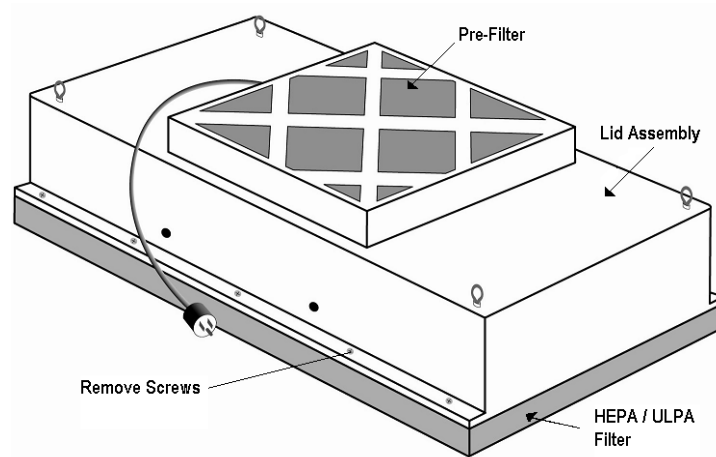



Removing the Fan/Filter Unit



Disconnect the unit from the electrical power source before attempting any service.

1. Disconnect the fan/filter unit from the power strip and roll up the power cord.
2. Attach double stick tape to all four sides of a polypropylene panel, as shown at right.
3. Place the polypropylene panel over the filter screen, making sure that it only adheres to the screen without overlapping the edges.
4. Push one side of the FFU up, rotate it 90° and lower it through the ceiling grid.
5. After replacing the filter, reverse these steps to reinstall the fan/filter unit in the ceiling grid.

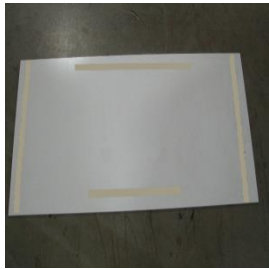


Figure 1



Figure 2



Figure 3



Figure 4

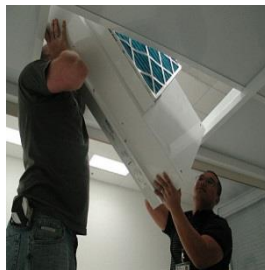


Figure 5



Figure 6



Cleanroom Maintenance

Useful Tools

Wipes

Wipes are used more frequently than any other cleaning product or tool. Selection of wipes should be based on intended usage. When selecting wipes you should consider things such as particle-shedding properties, chemical residue of the wiper content, static properties, absorbency and size. When using wipes, wipe in one direction from top to bottom or left to right. Use only slight overlapping strokes. Remove surface spots with commercial cleaner and woven polyester wipes. If cleaning with isopropyl alcohol or a similar cleaning agent, preform a full wipe-down of the sanitized area with deionized water.

Cavicide and Caviwipes and similar products contain ethylene glycol and benzethonium chloride, which are both caustic chemicals that will corrode stainless steel if not rinsed off with water afterwards. Any user disinfecting with Cavicide or similar product should preform a full wipe-down of the area with deionized water and dried.



WARNING: Always check chemical compatibility before cleaning plastic surfaces. Although vinyl and polyurethane withstand exposure to a wide range of common cleaning agents, repeated exposure to strong chemicals can cause damage.

Vacuums

There are a variety of different Vacuums available for your cleanroom. Selection of a vacuum will depend heavily on the application and the type of cleanroom you have. With all different types of sizes and filtration systems, select the one you feel would best suit the cleaning needs of your room. Refer to the Parts & Accessories section. For more information log on to our website at Terrauniversal.com

Mini-Environment Cleaning Kits

The ITW Texwipe Mini Environment Cleaning Kits are ideal for cleaning corners and difficult-to-reach locations inside the cleanroom. The kits include a cleaning tool (18" and 24" handles, 1 polyester foam pad, and 6 mop covers), one production bag of dry and pre-wetted wipers and an informational brochure with instructions on how to clean your equipment.

Designed to facilitate cleaning, the mop head has a low, flat profile with rounded corners and is autoclaveable. The swivel joint allows the user to reach inaccessible areas and replaceable foam pad ensures that the mop cover conforms to the surfaces that are being cleaned. The polyester knit fabrics used for the wipers and mop covers will not contaminate isolator surfaces when used in cleaning and disinfection operations.



5.0 Warranty

Products Manufactured by Terra: Terra Universal, Inc., warrants products that it manufactures to be free from defects for a period of 12 months for parts and 90 days for labor, commencing from the date of shipment. This limited warranty covers parts and labor, but not transportation and insurance charges. Terra's sole responsibility is to repair or replace, at its option, any part of the product that proves defective or malfunctioning during this time limit. In some cases, components incorporated in Terra Universal products are covered by additional warranties from component manufacturers; obtain specific information from Terra sales representatives. Repairs may be completed by 3rd party service agents approved by Terra Universal. Terra Universal reserves the rights to limit this warranty based on a service agent's travel, working hours, the site's entry restrictions and unobstructed access to serviceable components of the product. This warranty is void if the equipment is abused or modified by the customer, is operated outside Terra's operating instructions or specifications, or is used in any application other than that for which it is specified. This warranty does not include routine maintenance or service procedures, breakage of quartz baths after 60 days, shipping damage, nor damage from misuse, intentional or unintentional abuse, neglect, natural disasters, or acts of God.

Products Manufactured by Others: Terra Universal, Inc., warrants that, to the best of its ability, Terra's representations of products that are manufactured by others reflect the manufacturer's representations, subject to change without notice. Sole warranty for these products is the original manufacturer's warranty that is passed forward to the purchaser and constitutes the customer's sole remedy for these products. Detailed warranties for distributed products are available through Terra sales representatives.

Freight Shortage or Damage: Upon receipt of any equipment from Terra Universal, Inc., customer shall immediately unpack and inspect for damage or shortage. The customer shall not accept a damaged package or a short shipment until the carrier makes a "damage or shortage" notation on both the carrier's and customer's copy of the freight bill or delivery receipt. Service title passes when the shipment is loaded, so customer is responsible for filing and collecting a freight claim. Any replacement products must be ordered and paid for separately. For Terra's "Policy and Procedures for Returning Goods," see Terra's Internet site: www.TerraUniversal.com.

Generally, customers can improve the chance of collecting on a freight claim by following these procedures: 1) formally requesting that the carrier inspect the shipment immediately upon suspecting damage or shortage to verify condition; 2) notifying the carrier upon discovery of concealed damage and requesting an inspection within 15 days of receipt, both in person or phone and following up via mail; 3) keeping the shipment as intact as possible, including retaining original packaging materials and keeping the product as close to the original receiving location as possible; 4) holding salvage for disposition by the carrier.

All Claims: Terra Universal expressly disclaims all other warranties, expressed or implied or implied by statute, including the warranties of merchantability or fitness for intended use. Terra Universal is not responsible for consequential or incidental damages arising out of the purchase or use of the products supplied by Terra Universal. Terra Universal is not liable for damage to facilities, other equipment, products, property or personnel of others, or of their agents, suppliers, or affiliated parties, which is caused or alleged to have been caused by products supplied by Terra Universal. In any event or series of events, Terra Universal's total liability for any and all damages whatsoever is limited to the lesser of the actual damages or the original invoice cost of the items alleged to have caused the damage. The customer's sole and exclusive remedy for any cause of action whatsoever is repair or replacement of the non-conforming products or refund of the actual purchase price, at the sole option of Terra Universal. All claims must be made in writing within 90 days of the date the product was shipped. Any claims not made within this time limit shall be deemed waived by the customer. Terra Universal is not responsible for any additional costs of repair caused by poor packaging or in-shipment damage during return.

Warranty Returns: All warranty returns must be authorized in advance by Terra Universal and approved under an RMA. Unless approved in advance for good reason, all returns must be in original condition, including all manuals, and must be packaged in original packaging materials. All returned goods are to be shipped to Terra Universal, freight prepaid at customer's expense. See Terra's "Policy and Procedure for Returned Goods."

Terra Universal Warranty Policy: <https://www.terrauniversal.com/warranty/>

*Thank you for ordering from
Terra Universal!*



6.0 Spare Parts & Accessories

For replacement parts and accessories for your cleanroom order at
Terra Universal.com or Call 714-578-6000

Filter Replacement

Replacement Filters for Terra WhisperFlow™ Fan Filter Units		
Type	Size	TUI Part #
HEPA	2 X 2	6601-27
	2 X 3	6601-26
	2 X 4	6601-25
ULPA	2 X 2	6601-30
	2 X 3	6601-29
	2 X 4	6601-28

Cleanroom and Laboratory Wipers



Type	Size	Qty	TUI Part #
Cotton Wiper	9" x 9"	150/bag	5605-07
Cotton Wiper	12" x 12"	75/bag	5605-02

White twill-patterned wiper, made of 100% long-staple cotton is tightly woven for strength and durability, while also clean and absorbent. Fabric is bias-cut and cross section of 120x60 threads per square inch. The woven twill pattern makes this wiper ideal for cleaning, polishing and burnishing magnetic disk media surfaces. Ideal in high heat applications, and absorbing aqueous and organic solvent spills. Suitable for Class 1000 and higher cleanrooms.

Laundered Polyester Wipers



Size	Standard 209E	Qty	TUI Part #
9" x 9"	Class 1	150/bag	5605-00
12" x 12"	Class 10	75/bag	5605-08

Laundered wipers are ideal for critical environments where ultra-low particle/fiber generation and absorbent capacity is important. Soft yet durable construction cleans surfaces without scratching.



Softwall Modular Cleanroom

Caster



Type	TUI Part #
6" Locking Caster	6704-08

This 6 " locking stainless steel swivel casters feature nonshedding polyurethane wheels. Capacity: 300 lbs. each. Order one per vertical post. Casters raise the frame 7.125". If casters are ordered, the frame will be made shorter to keep the floor to filter face dimension to specification.

MicroVac Portable Cleanroom Vacuum Cleaner



Type	TUI Part #
110 Vac/60Hz	5100-00
220 Vac/50Hz	5100-00-220

The MicroVac™ is a portable vacuum cleaner suitable for use in cleanroom environments. It incorporates a four-stage filter system to remove particles and dust from the air before returning the air to the workplace. The small size, light weight, and flexible attachments of this cleaner make it ideal for cleaning tight corners and delicate equipment that a larger cleaner just can't reach!

HEPA-Filtered Vacuum Cleaner



Type	TUI Part #
110 Vac/60Hz	1001-00

This HEPA-Filtered Vacuum Cleaner combines the power and capacity for heavy-duty operation with a lightweight, compact design that makes it as convenient as it is effective. The unit's HEPA (High Efficiency Particulate Air) filter removes 99.999% of all particles 0.3µm or larger, exceeding federal standards and accepted HEPA filtration levels.

ULPA Filtration Vacuum Cleaner



Type	TUI Part #
115 Vac/60Hz	1764-00
220 Vac/50Hz	1764-00-220

This Cleanroom Vacuum Cleaner is used for cleaning the workplace or picking up spills of hazardous dust and particles. Its unique four-stage filter system removes particles 0.12 microns in size at 99.9995% efficiency, before returning the air to the workplace.



Softwall Modular Cleanroom

Light Panel



Type	TUI Part #
120-277V Vac. 50/60 Hz	3800-41B

Yellow 12-foot Power Cables

For connection of light or fan filter unit to Power Distribution Module (electrical box)



No. EL01298: for Fan Filter Units
4 Pin – up to 600 VAC



No. EL01297: for Light Panels
3 Pin – Up to 600 VAC